# Manual Rotor Beater Mill SR 300









# Copyright

© Copyright by Retsch GmbH Retsch-Allee 1-5 42781 Haan Germany



# **Table of Contents**

1		Note	es on the Manual	. 6
	1.1		Disclaimer	. 6
	1.2		Copyright	
	1.3		Explanations of the Safety Instructions	
	1.4		General Safety Instructions	
	1.5		Repairs	
2			firmation Form for the Managing Operator	
3			nnical Data	
	3.1		Protective Equipment	
	3.2		Degree of Protection	
	3.4		Electromagnetic Compatibility (EMC)	
	3.5		Rated Power	
	3.6		Motor Rotation Speed	
	3.7		Dimensions and Weight	
	3.8		Required Floor Space	
	3.9 3.10		Receptacle Volume	
	3.11		Feed Grain SizeSuitable Sieve Diameters	
	-			
4			kaging, Transport and Installation	
	4.1		Packaging	
	4.2 4.3		ransport  Temperature Fluctuations and Condensation	
	4.3		Conditions for the Installation Site	
	4.5		Electrical Connection	
	4.6		Type Plate Description	
5		Firef	t Commissioning	18
•	5.1		Setting up as Table-top Device	
	5.2		Setting up on Base Frame	
c			rating the Device	
6		•		
	6.1 6.2		Use of the Device for the Intended Purpose	
	6.3		/iews of the Instrument	
		.3.1	Front	
	-	.3.2	Back	
	6.4		Switching On / Off	
	6.5		Opening and Closing of the Device	
	6.	.5.1	Opening	
	6.	.5.2	Closing	
	6.6	Е	Emergency Unlocking	31
	6.7	lı	nserting the Grinding Set	32
	6.	7.1	Inserting the Grinding Chamber Cartridge	33
		7.2	Inserting the Rotor	
		.7.3	Inserting the Retaining Frame	
	6.8		Removing the Grinding Set	
	6.9	Λ	Mounting the Sample Receptacle	35
7		Con	trolling the Device	37
	7.1	C	Operating Controls, Display and Functions	37
	7.2		Nanual Mode	
	7.	.2.1	Setting the Rotation Speed	
		2.2	Start Process	
	7.	.2.3	Feeding Sample Material	38



7.2.4 Stop Process	38
8 Error Messages and Information Notes	39
8.1 Error Messages	39
9 Return for Service and Maintenance	
10 Cleaning, Wear and Maintenance	
10.1 Cleaning	41
10.1.1 Cleaning the Grinding Set	42
10.1.2 Cleaning the Feed Hopper	
10.2 Wear	44
10.3 Maintenance	
11 Accessories	46
11.1 Vibratory Feeder DR 100	46
11.2 Cyclone Separator	48
12 Disposal	54
13 Index	55





## 1 Notes on the Manual

#### Dear user,

please read the following manual referring to this device carefully before starting any installation, commissioning and operation.

This manual is a technical guide on how to operate the device safely and it contains all the information required for the areas specified in the table of contents. This technical documentation is a reference and instruction manual. The individual chapters are complete in themselves. Familiarity (of the respective target groups defined per area) with the relevant chapters is a precondition for the safe and appropriate use of the device.

This manual does not contain any repair instructions. In case of any obscurities or questions with regards to this document or the device, as well as if faults arise or repairs are necessary, please contact your supplier or get in touch with Retsch GmbH directly.

Application-technical information relating to samples to be processed are not or only to a certain extend included. However, more information thereof can be found in the internet on the webpage of the respective device on the Retsch GmbH homepage (http://www.retsch.com).

#### **Revision status:**

This document revision 0002 refers to the manual "Rotor Beater Mill SR 300" in compliance with the Directive of Machinery 2006/42/EC.

#### 1.1 Disclaimer

This document has been prepared with due care. Technical and software based modifications are reserved. No liability is assumed for data loss, personnel injury or damage to the device which results from the failure to observe the instructions and/or warnings in this document.

# 1.2 Copyright

This document or parts of it or its content may not be reproduced, distributed, edited or copied in any form without prior written permission of Retsch GmbH. Damage claims shall be asserted in the case of infringements.



# 1.3 Explanations of the Safety Instructions

In this document the following signs and symbols are being used:

<ul><li>①</li><li>→</li></ul>	Reference to a recommendation and/or an important information		
$\rightarrow$	Reference to a chapter, table or figure		
$\Rightarrow$	Action instruction		
Name	Software menu function		
[Name]	Software button		
(Name)	Software checkbox		

In this document the following safety instructions warn of possible dangers and damages:



D1.0000

### Type of danger / personal injury

Source of danger

- Possible consequences if the dangers are not observed.
- · Instructions and information on how the dangers are to be avoided.

**Life-threatening personal injuries** may result from disrespecting the safety instruction for danger. There exists a very high risk of hazard of life or permanent injury to personnel. Additionally, in continuous text or action instructions the signal word **A DANGER** is used.



#### WARNING

W1.0000

## Type of danger / personal injury

Source of danger

- Possible consequences if the dangers are not observed.
- . Instructions and information on how the dangers are to be avoided.

**Serious personal injuries** may result from disrespecting the warning notice. There exists an elevated risk of an accident or severe injury to personnel. Additionally, in continuous text or action instructions the signal word **A WARNING** is used.



#### CAUTION

C1.0000

### Type of danger / personal injury

Source of danger

- Possible consequences if the dangers are not observed.
- Instructions and information on how the dangers are to be avoided.

**Moderate or mild personal injuries** may result from disrespecting the safety instruction for caution. There exists a medium or low risk of an accident or injury to personnel. Additionally, in continuous text or action instructions the signal word **A CAUTION** is used.



**NOTICE** 

N1.0000

#### Type of property damage

Source of property damage

- Possible consequences if the notices are not observed.
- Instructions and information on how the property damages are to be avoided.

**Property damages** may result from disrespecting the notice. However, there exists no risk of an injury to personnel. Additionally, in continuous text or action instructions the signal word **NOTICE** is used.

# 1.4 General Safety Instructions



#### **CAUTION**

C2.0002

#### Read the manual

Non-observance of the operating instructions

- The non-observance of this manual can result in personal injuries.
- · Read the manual before using the device.



#### Target group:

All persons concerned with this device in any form.

This device is a modern, high performance product from Retsch GmbH and complies with the state of the art. Operational safety is given if the device is handled for the intended purpose and attention is given to this technical documentation.

#### Safety manager:

The managing operator himself must ensure that the people entrusted with working on the device...

- have noted and understood all the regulations regarding safety,
- are familiar before starting work with all the operating instructions and specifications for the target group relevant to their work,
- have unrestricted and free access to the technical documentation of this device,
- are familiarising new personnel with the safe handling of the device and its use for its intended purpose before starting work either by verbal instructions from a competent person and/or by means of this technical documentation.

▲ CAUTION Improper operation can result in personal injuries and material damage. The managing operator himself is responsible for his own safety and that of his employees. The managing operator himself is responsible that no unauthorised person has access to the device.

▲ CAUTION Persons who are under the influence of intoxicating substances (pharmaceuticals, drugs, alcohol), fatigue or health disorders must not operate the device.





# **CAUTION**

C3.0015

#### Changes to the device

Improper modifications

- Changes to the device can lead to personal injuries.
- Do not make any modification to the device.
- Use spare parts and accessories that have been approved by Retsch GmbH exclusively.

## **NOTICE**

N2.0012

#### Changes to the device

Improper modifications

- The conformity declared by Retsch GmbH with the European Directives will lose its validity.
- Any warranty claims will be terminated.
- Do not make any modification to the device.
- Use spare parts and accessories that have been approved by Retsch GmbH exclusively.

# 1.5 Repairs

This manual does not contain any repair instructions. For safety reasons, repairs may only be carried out by Retsch GmbH or an authorised representative or by qualified service technicians.

#### In case of repair, please inform...

- ...the Retsch GmbH representative in your country,
- ...your supplier, or
- ...Retsch GmbH directly.

Service address:					



# **2** Confirmation Form for the Managing Operator

This manual contains essential instructions for operating and maintaining the device which must be strictly observed. It is essential that they be read by the user and by the qualified staff responsible for the device before the device is commissioned. This manual must be available and accessible at the place of use at all times.

The user of the device herewith confirms to the managing operator (owner) that he has received sufficient instructions about the operation and maintenance of the system. The user has received the manual, has read and taken note of its contents and consequently has all the information required for safe operation and is sufficiently familiar with the device.

The managing operator should for legal protection have the user confirm the instruction about the operation of the device.

I have read and taken note of the contents of all chapters in this manual as well as all safety instructions and warnings.	
User	
Surname, first name (block letters)	
Position in the company	
Place, date and signature	
Managing operator or service technician	
Surname, first name (block letters)	
Position in the company	
Place, date and signature	



#### 3 Technical Data

# 3.1 Protective Equipment

- This device is equipped with a door lock. The locking mechanism prevents the device from being started when in an unsafe state.
- The device can only be started when the door is closed.
- It is only possible to open the door when the device has come to a halt.

## 3.2 Degree of Protection

- IP41

#### 3.3 Emissions



C4.0020

#### Possibility of acoustic signals not being heard

Loud grinding noises

- Possible acoustic alarms and voice communication might not be heard.
- Consider the volume of the grinding noise in relation to other acoustic signals in the work environment. Additional visual signals may be used.

# **A** CAUTION

C5.0044

#### **Hearing damage**

A high sound level may be generated depending on the type of material, the grinding set used, the grinding frequency set and the duration of the grinding



- Excessive noise in terms of level and duration can cause impairments or permanent damage to hearing.
- Ensure that suitable noise protection measures are taken, or wear ear protection.

#### Sound parameters:

The sound parameters are influenced by the properties of the sample material and the rotation speed.

#### Example 1:

Feed material:	Limestone (< 5 mm)
Bottom sieve:	0.5 mm
Speed:	3 000 rpm

At these operating conditions, the workplace related equivalent continuous sound level  $L_{eq} = 76.3 \text{ dB(A)}$ .



#### Example 2:

Feed material:	Limestone (< 5 mm)
Bottom sieve:	0.5 mm
Speed:	5 000 rpm

At these operating conditions, the workplace related equivalent continuous sound level  $L_{eq} = 72.8 \text{ dB(A)}$ .

#### Example 3:

Feed material:	Limestone (< 5 mm)
Bottom sieve:	0.5 mm
Speed:	10 000 rpm

At these operating conditions, the workplace related equivalent continuous sound level  $L_{eq} = 77.2 \text{ dB(A)}$ .

# 3.4 Electromagnetic Compatibility (EMC)

- EMC class according to DIN EN 55011: B

#### 3.5 Rated Power

~ 2500 W (VA)

# 3.6 Motor Rotation Speed

- Rated motor rotation speed: 3 000 10 000 revolutions per minute (rpm).
- Adjustable in increments of 500 rpm.

In order to reduce the heating of the grinding material caused by high friction forces at high speeds, a distance rotor can be used for the grinding of heat-sensitive products.

# 3.7 Dimensions and Weight

All specifications, unless otherwise stated, include the optional base frame.

Height: 1 200 mm
Height incl. vibratory feeder DR 100: ~ 1 450 mm
Width: 580 mm
Width incl. cyclone separator: 930 mm
Depth: 700 mm
Weight: ~ 60 kg
Weight without base frame: ~ 50 kg



## 3.8 Required Floor Space

# **A** CAUTION

C6.0007

#### Falling down of the device

Incorrect positioning or insufficient working space

- Due to its weight, the device can inflict personal injury if it falls down.
- The device must, when used as table-top device, only be operated on a sufficiently large, strong and stable workplace.
- The device must, when used as table-top device, be firmly screwed to the working surface.

Width with open door: ~ 740 mmDepth with open door: ~ 640 mm

Width of the base: 940 mm (without cyclone separator)

1 130 mm (with cyclone separator)

Depth of the base: 740 mm (without base frame)

800 mm (with base frame)

#### **Location requirements:**

When used as table-top device, the device must be bolted to a vibration-free, plane, stable and free surface. When used as floor-standing device (mounted on the optional base frame), the device must be placed on a level base and secured against rolling away.

# 3.9 Receptacle Volume

The receptacle volume (feed volume) depends on the sample material and the collecting receptacle used.

Feed volume: < 5 I (with 5 I collecting receptacle)</li>

max. 26 I (with 30 I collecting receptacle)

#### 3.10 Feed Grain Size

Feed grain size: ≤ 25 mm

## 3.11 Suitable Sieve Diameters

Trapezoidal perforation: 0.08 mm / 0.12 mm / 0.20 mm / 0.25 mm / 0.35 mm / 0.50 mm /

0.75 mm / 1.00 mm / 1.25 mm / 1.50 mm / 2.00 mm

Round perforation: 3.00 mm / 4.00 mm / 10.00 mm

All sieve inserts are available as 360° sieve insert (ring sieve), as well as 180° sieve insert. The 360° sieve inserts are particularly suitable for the grinding of soft and brittle materials. To achieve an additional size reduction effect through impact, a 180° grinding insert can be used together with a 180° sieve insert for medium-hard materials.



# 4 Packaging, Transport and Installation

# 4.1 Packaging

The packaging has been adapted to the mode of transport. It complies with the generally applicable packaging guidelines.

### **NOTICE**

N3.0001

#### Storage of the packaging

- In the event of a complaint or return, the warranty claim may be endangered if the packaging is inadequate or the device has not been secured correctly.
- Keep the packaging for the duration of the warranty period.

# 4.2 Transport



#### **WARNING**

W2.0005

#### Serious personal injury

#### Falling loads



- Due to the heavy weight of the device, serious personal injuries can be caused if it falls down.
- Lifting above head height is not permissible!

#### **NOTICE**

N4.0017

#### **Transport**

- Mechanical or electronic components may be damaged.
- The device may not be knocked, shaken or thrown during transport.

#### **NOTICE**

N5.0014

#### **Complaints**

Incomplete delivery or transportation damage

- The forwarding agent and Retsch GmbH must be notified immediately in the event of transportation damage. It is otherwise possible that subsequent complaints will not be recognised.
- Please check the delivery on receipt of the device for its completeness and intactness.
- Notify your forwarding agent and Retsch GmbH within 24 hours.





Fig. 1: Transportation lug

▲ CAUTION The weight of the SR 300 without collecting receptacle and base frame amounts approx. 50 kg. The device may only be lifted or transported by at least two people.

The SR 300 has a transportation lug (**A**) at which the device can be lifted and transported by means of hoists.

- ⇒ Attach the hoist only to the supplied transportation lug (**A**).
- ⇒ Please insert a piece of cloth, polystyrene or cardboard between the rear edge of the feed hopper and the sling or hoisting chain, in order to avoid scratches on the feed hopper.
- ⇒ Lift the device slowly and stabilise it to prevent tipping backwards.

**A** CAUTION Only use suitable hoist that is designed for the weight of the device.

If desired, the transportation lug can be removed after installation and the threaded hole can be sealed with the supplied plastic screw (**PS**).

#### 4.3 Temperature Fluctuations and Condensation

## **NOTICE**

N6.0016

#### **Temperature fluctuations**

The device may be subject to strong temperature fluctuations during transport (e.g. aircraft transport)

- The resultant condensed water may damage electronic components.
- Wait before commissioning until the device has been acclimatised.

#### Temporary storage:

Also in case of an interim storage the device must be stored dry and within the specified ambient temperature range.

#### 4.4 Conditions for the Installation Site

Installation height: max. 2 000 m above sea level

Ambient temperature: 5 °C – 40 °C



**NOTICE** 

N7.0021

#### Ambient temperature

Temperatures outside the permitted range

- Electronic and mechanical components may be damaged.
- The performance data alter to an unknown extent.
- Do not exceed or fall below the permitted temperature range (5 °C to 40 °C ambient temperature) of the device.
- Maximum relative humidity < 80 % (at ambient temperatures ≤ 31 °C)</li>

For ambient temperatures  $U_T$  between 31 °C and 40 °C, the maximum relative humidity value  $L_F$  linearly decreases according to  $L_F = -(U_T - 55) / 0.3$ :

Ambient temperature	Max. rel. humidity
≤ 31 °C	80 %
33 °C	73.3 %
35 °C	66.7 %
37 °C	60 %
39 °C	53.3 %
40 °C	50 %

**NOTICE** 

N8.0015

#### Humidity

High relative humidity

- Electronic and mechanical components may be damaged.
- The performance data alter to an unknown extent.
- The relative humidity in the vicinity of the device should be kept as low as possible.

#### 4.5 Electrical Connection

**WARNING** When connecting the power cable to the mains supply, use an external fuse that complies with the regulations applicable to the place of installation.

- Check the type plate for details on the necessary voltage and frequency for the device.
- The listed values must agree with the existing mains supply.
- Only use the supplied power cable to connect the device to the mains supply.

**NOTICE** 

N9.0022

#### **Electrical connection**

Failure to observe the values on the type plate

- Electronic and mechanical components may be damaged.
- Connect the device only to mains supply matching the values on the type plate.



# 4.6 Type Plate Description

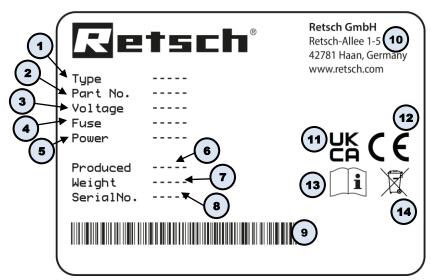


Fig. 2: Type plate

- 1 Device designation
- 2 Part number
- 3 Power version, Mains frequency
- 4 Fuse type and fuse strength
- 5 Capacity, Amperage
- 6 Year of production
- 7 Weight
- 8 Serial number
- 9 Bar code
- 10 Manufacturer's address
- 11 UKCA marking
- 12 CE marking
- 13 Safety warning: Read the manual
- 14 Disposal label
- ① In the case of queries please provide the device designation (1) or part number (2), as well as the serial number (8) of the device.



# 5 First Commissioning

# A

#### **WARNING**

W3.0002

#### Danger to life through electric shock

Damaged power cable

- An electric shock can cause burns, cardiac arrhythmia, respiratory arrest, as well as cardiac arrest.
- Never use a damaged power cable to connect the device to the mains!
- Check the power cable and the plug for any damage before use.



## **NOTICE**

N10.0002

#### Setting up the device

Disconnecting the device from the mains

- A separation of the device from the mains must be possible at any time.
- Set up the device in such a way, that the connection for the power cable is always easily accessible.

## **NOTICE**

N11.0004

#### Setting up the device

Vibrations during operation

- Depending on the operating mode of the device, slight vibrations may occur.
- Set up the device only on a vibration-free, plane and stable surface.



# 5.1 Setting up as Table-top Device

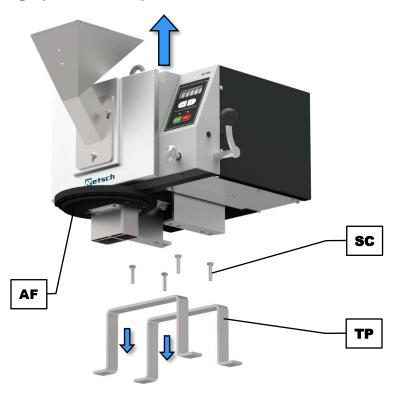


Fig. 3: Removing the transport profiles

- ⇒ Loosen the four hexagon socket head screws (**SC**) to remove the transport profiles (**TP**) from the SR 300.
- ⇒ Position the SR 300 on the designated table as close as possible to the front edge of the table. Make sure that the discharge flange (**AF**) for the sample receptacle does not touch the edge of the table.

**NOTICE** The table must be designed for the weight of approx. 50 kg of the SR 300! Furthermore, the table must be plane, stable and vibration free!



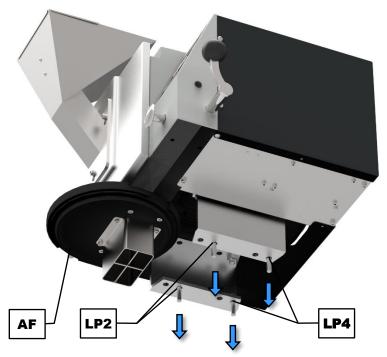


Fig. 4: Table mounting

⇒ Bolt the SR 300 to the tabletop by means of four M8 screws of the required length (not supplied). Use the second (LP2) and fourth pair of holes (LP4) of the U-profiles for this purpose.

# 5.2 Setting up on Base Frame

① The base frame for the SR 300 is available as an optional accessory from your supplier or from Retsch GmbH directly.

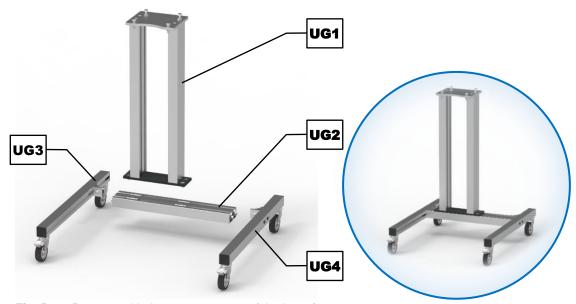


Fig. 5: Pre-assembled component parts of the base frame

The following component parts are already pre-assembled for an easier assembly of the base frame:



- base (**UG1**)
- cross bar (UG2)
- left roller rail (UG3)
- right roller rail (UG4)

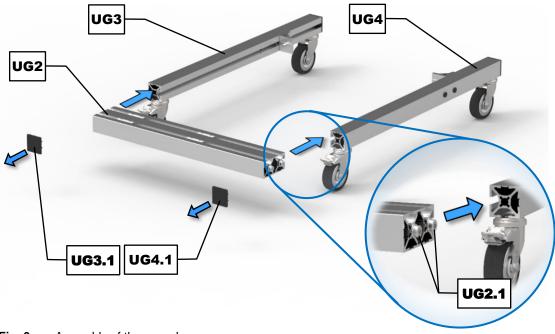


Fig. 6: Assembly of the cross bar

- ⇒ Remove the two front covers (**UG3.1**) and (**UG4.1**) from the left and right roller rails (**UG3**) and (**UG4**).
- ⇒ Loosen the four hexagon socket head screws (**UG2.1**) on the cross bar (**UG2**) by means of the supplied hexagon socket wrench (**IM**).
- ⇒ Slide the cross bar (**UG2**) into the lateral guides of the left and right roller rails (**UG3**) and (**UG4**). When doing so, make sure that the washers of the four hexagon socket head screws (**UG2.1**) are located inside the roller rails.

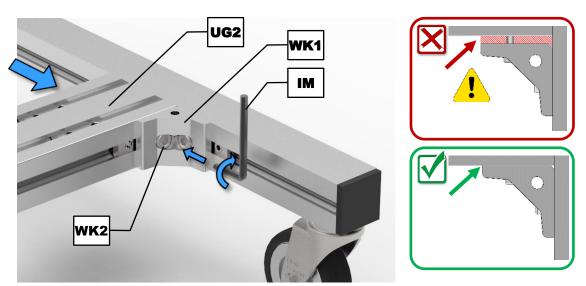


Fig. 7: Attaching the cross bar to the brackets



- ⇒ Use the two supplied M8x16 hexagon socket head screws (**WK2**) to screw the cross bar (**UG2**) to the two brackets (**WK1**) of the left and right roller rails.
- ⇒ Tighten the hexagon socket head screws (WK2) firmly to the left and right brackets (WK1).

**NOTICE** Ensure that the cross bar (**UG2**) lies flush against the two pre-assembled brackets (**WK1**).

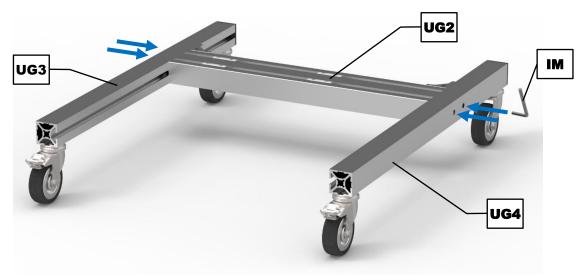


Fig. 8: Screwing tight the cross bar

⇒ Tighten again the four hexagon socket head screws (**UG2.1**) of the cross bar (**UG2**) firmly from the side of both roller rails.

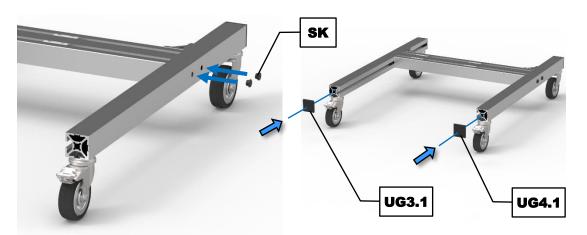


Fig. 9: Installing the protective caps and covers

- ⇒ Place the four supplied black protective caps (**SK**) on the lateral openings of the two roller rails
- ⇒ Reinstall the two front covers (UG3.1) and (UG4.1) of the left and right roller rail.



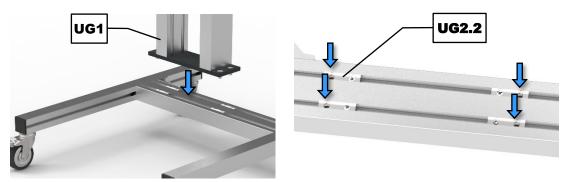


Fig. 10: Positioning the base

⇒ Place the base (**UG1**) on the cross bar (**UG2**) and use the four supplied M8x20 hexagon socket countersunk head screws (**UG1.1**) to tighten the base.

**NOTICE** The base should be positioned in such a way, that the left edge of the base is located 90 mm from the outer edge of the left roller rail (**UG3**).

⇒ If necessary, adjust the position of the sliding blocks (UG2.2).

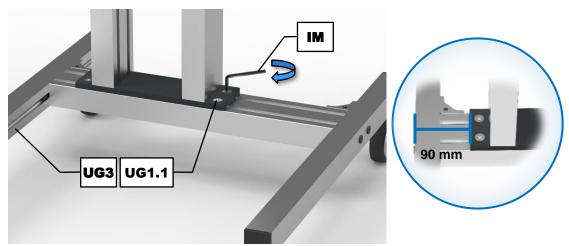


Fig. 11: Screwing tight the base

- ⇒ Place the SR 300 on the base frame.
- ⇒ Use the four supplied M8x35 hexagon socket head screws (**UG1.2**), including the locknuts and washers, to screw the U-profiles of the SR 300 tight on the base (**UG1**).



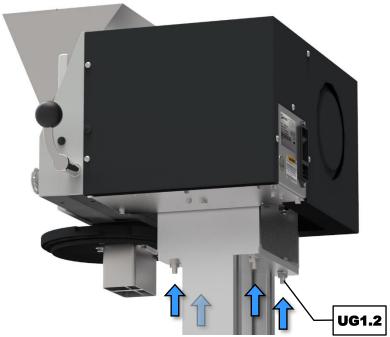


Fig. 12: Mounting the SR 300



# 6 Operating the Device

## 6.1 Use of the Device for the Intended Purpose

# A

#### **CAUTION**

C7.0005

#### Risk of explosion or fire

Potentially explosive atmosphere

- On account of its design, the device is not suitable for use in potentially explosive atmospheres.
- Do not operate the device in a potentially explosive atmosphere.

# A

#### CAUTION

C8.0006

#### Danger of personal injury

Hazardous sample material

- Depending on the dangerous nature of the sample material necessary measures must be taken to rule out any danger of personal injury.
- Observe the material safety data sheets of the sample material.



# A

#### **CAUTION**

C9.0004

#### Risk of explosion or fire

Changing sample properties



- The properties and therefore also the hazardousness of the sample can change during the grinding process.
- Do not use any substances in this device which carry the risk of explosion or fire.
- Observe the material safety data sheets of the sample material.


# A

#### **CAUTION**

C10.0010

#### **Chemical reactivity**

Changing sample properties

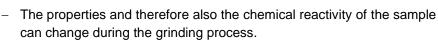

- Do not use any substances in this device which carry the risk of increasing chemical reactivity during the grinding process.
- Observe the material safety data sheets of the grinding material.



This Rotor Beater Mill of the Retsch GmbH is a laboratory device. It is used for the batch-wise or continuous coarse and fine grinding of dry, soft to medium-hard materials such as soils, bentonite, chemical products, fertilizers, drugs, feed pellets, spices, gypsum, graphite, resins, coal, coke, herbs, plastics, food, plant parts, pharmaceutical products, powder lacquer, rice, seeds, salts, sugar and many other substances. The device allows for a particularly effective grinding of heterogeneous material mixtures. The analytical fineness is usually achieved in a single step. Thereby, the grinding material is hardly heated, so that the SR 300 is also suitable for temperature-sensitive materials. Depending on the mesh size of the sieve and the fracture properties of the grinding material, the achievable final fineness can be < 50  $\mu$ m.



The Rotor Beater Mill of the Retsch GmbH is universally deployed in almost all areas of industry and research, from sample preparation in the laboratory to larger sample volumes in a technical centre or plant.

Only grinding sets from Retsch GmbH may be used. This device is generally not designed for the grinding of wet or moist materials.

## **NOTICE**

N12.0007

## Range of application of the device

Long-term operation

- This laboratory device is designed for eight-hour single-shift operation with a duty cycle of 30 %.
- This device may not be used as a production machine nor is it intended for continuous operation.

# 6.2 Principle of Operation

The grinding and deagglomeration in the SR 300 is achieved by impact, collision and shearing effects. The feed material passes through the feed hopper into the grinding chamber where it is comminuted between the rotor and the selected ring sieve or sieve and grinding insert. As soon as the grinding material is smaller than the mesh size of the sieve, it enters the collecting receptacle.



# 6.3 Views of the Instrument

# 6.3.1 Front

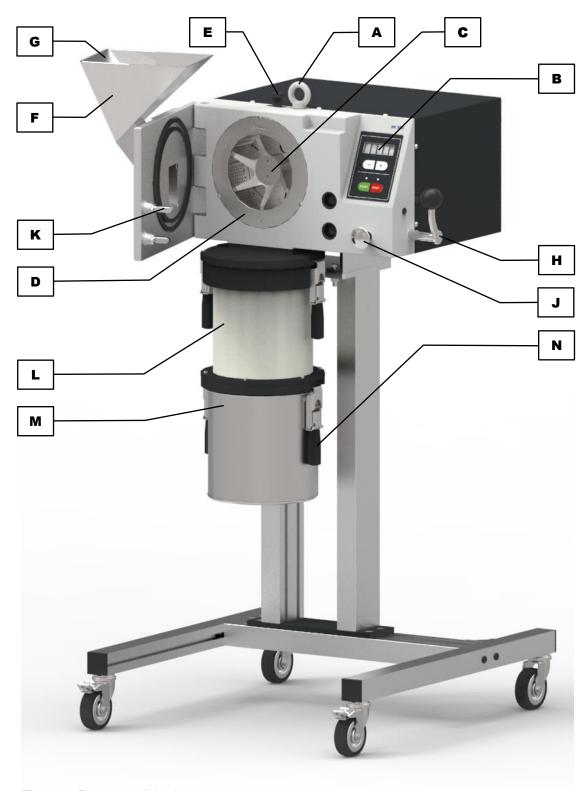


Fig. 13: Front view of the device



Element	Description	Function
A	Transportation lug	Attachment point for hoist
В	Operating controls	Operation of the device
С	Rotor	Crushes the sample material
D	Retaining frame with sieve insert	Sieves the grinding material
E	Locking pin	Holds the grinding chamber cartridge in position
F	Feed hopper	Takes the sample material
G	Splashback protection	Prevents material ejection
Н	Locking lever	Locks or opens the door
J	Catch mechanism	Locks the locking lever
K	Locking mechanism	Keeps the device closed
L	Filter hose	Ensures pressure compensation
M	Collecting receptacle	Receives the crushed grinding material
N	Clamping lever	Seals the collecting receptacle or the filter hose. Also serves to carry the collecting receptacle in unfolded state

# 6.3.2 Back



Fig. 14: Back view of the device



Element	Description	Function
Α	Transportation lug	Attachment point for hoist
E	Locking pin	Holds the grinding chamber cartridge in position
F	Feed hopper	Takes the sample material
Н	Locking lever	Locks or opens the door
I	Mains switch	Switches the device on and off, disconnects the device from the mains
0	Type plate	Lists, among others, the voltage type, the serial number and the type of the device
P	Sticker "Manual"	Reminds to read the manual
R	Warning sign "Disconnect from the mains"	Warning of electric shock
S	Mains connection	Connection for the power cable

# 6.4 Switching On / Off

⇒ Turn on the SR 300 with the mains switch (I) on the back side of the device.

When the device is switched off, it is completely disconnected from the mains.

# 6.5 Opening and Closing of the Device

The SR 300 has a double locking mechanism (**K**) that locks the door both, mechanically and electrically. The electrical locking mechanism is closed by pressing the button, and by pressing the button opened again after the motor has come to a complete halt. The mechanical locking mechanism is closed or opened by actuating the locking lever (**H**) and the catch mechanism (**J**).



## 6.5.1 Opening



Fig. 15: Opening the device

- ⇒ Connect the device to the mains.
- ⇒ Switch the device on by means of the mains switch (I).

When the door is closed, the information note "H42" appears in the display. This informs the user to once open and close the door for safety reasons.

- ⇒ Pull and hold the catch mechanism (**J**) forwards.
- ⇒ Turn the locking lever (**H**) forwards. The locking mechanism (**K**) is released and the door can now be opened.

**NOTICE** If an attempt is made to open the device mechanically during operation, the error "E51" appears in the display. This indicates a fault in the safety switch.

- ⇒ Switch off the device by means of the mains switch (I).
- ⇒ Press the mains switch (I) again to turn the device on. "H42" appears in the display.
- ⇒ Open the door.

# 6.5.2 Closing

- ⇒ Close the door.
- ⇒ Turn the locking lever (**H**) backwards. The catch mechanism (**J**) snaps back into place by itself

If the door is not correctly locked, the information note "H41" appears in the display when the button is pressed. This instructs the user to close the door.



# 6.6 Emergency Unlocking

# **A** CAUTION

#### **Emergency unlocking**

Drive continuing to run

- Substantial risk of injury due to the drive and associated device parts running on a long time without being braked!
- Activate the emergency unlocking only when the device has come to a complete stop and when disconnected from the mains supply.

C11.0009

The emergency unlocking is used for manual unlocking of the electrical locking mechanism in the event of a power failure. To open the door, the mechanical locking mechanism must also be opened by pulling the catch mechanism (**J**) and turning the locking lever (**H**).

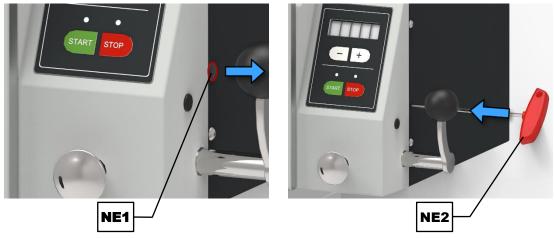


Fig. 16: Position of the emergency unlocking

- ⇒ Remove the protective cap (**NE1**) on the right side of the device.
- ⇒ Insert the supplied hexagon socket wrench (NE2) into the lateral opening.

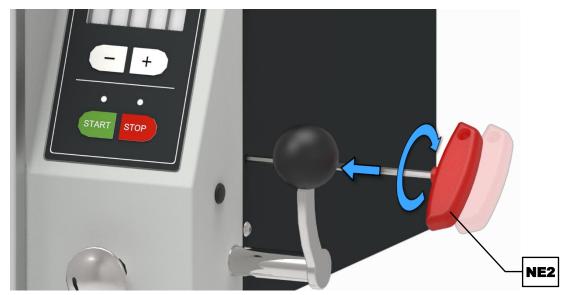


Fig. 17: Emergency unlocking of the device



To unlock the locking mechanism (**K**) the hexagon socket wrench (**NE2**) must be pressed further into the opening with a little force.

⇒ While pressing, turn the hexagon socket wrench (**NE2**) clockwise up to the stop.

The upper, electrical locking mechanism of the door is now unlocked and the door can be opened, if previously the lower, mechanical locking mechanism has been unlocked.

# 6.7 Inserting the Grinding Set



# Cutting injuries

Sharp cutting edges

- The sharp cutting edges on the rotors and/or the cutting bars in the grinding chamber can lead to hand lacerations.
- Grasp the rotor only on the hub and do not touch the cutting edges or bars.
- · Wear cut-resistant safety gloves.



C12.0085


N13.0066

# **NOTICE**

#### Wear or damage of the device

Operation without grinding set

- During operation of the device without grinding set, excessive wear or damage to the device may occur.
- Operate the device only with a grinding set mounted.

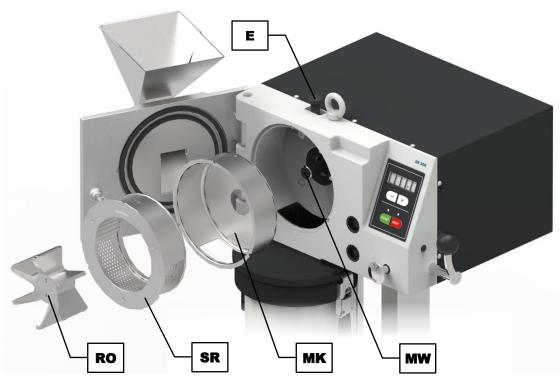

**Fig. 18:** Insertion of the grinding set consisting of grinding chamber cartridge, retaining frame and rotor



## 6.7.1 Inserting the Grinding Chamber Cartridge



Fig. 19: Inserting the grinding chamber cartridge

- ⇒ Pull up and hold the locking pin (**E**).
- ⇒ Align the grinding chamber cartridge so that the opening (MK1) for the locking pin is at the top, and the three recesses (MK2) are aligned with the corresponding cylinder pins.
- ⇒ Slide the grinding chamber cartridge into the grinding chamber.
- ⇒ Release the locking pin (**E**) to lock it into the opening (**MK1**) of the grinding chamber cartridge.

# 6.7.2 Inserting the Rotor



Fig. 20: Inserting the rotor

- ⇒ Use the supplied holding key (**HS**) to hold the motor shaft (**MW**) in place.
- ⇒ Grasp the rotor (RO) so that the rear hub (RO1) is aligned properly with the feather key (PF) on the motor shaft (MW).



⇒ Slide the rotor onto the motor shaft up to the stop.

**NOTICE** If it is hard to slide the rotor or if it cannot be pushed up to the stop, check the correct and firm fit of the feather key (**PF**) on the motor shaft (**MW**). In addition, the motor shaft can be oiled with some machine oil.

#### 6.7.3 Inserting the Retaining Frame

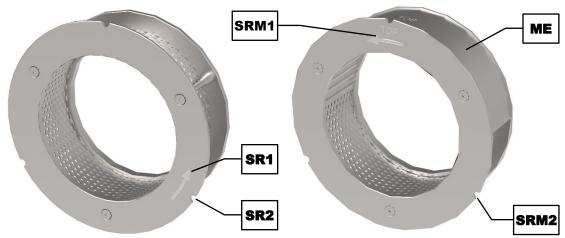


Fig. 21: Retaining frame for 360° sieve inserts (left) and 180° sieve inserts (right)

Both, the retaining frame for the 360° sieve inserts and the 180° sieve inserts have a direction arrow (**SR1**) or (**SRM1**) on the front, which indicates the direction of rotation of the rotor (counterclockwise).

- ⇒ Pull up and hold the locking pin (**E**).
- ⇒ Align the retaining frame so that the direction arrow (SR1) or (SRM1) is facing forward and the three recesses (SR2), (SRM2) are aligned with the corresponding cylinder pins. In the case of the retaining frame for the 180° sieve inserts, "TOP" must additionally be on top.
- ⇒ Slide the retaining frame into the grinding chamber cartridge (**MK**).
- ⇒ Release the locking pin (**E**).

#### Changing the sieve insert:

All available 360° and 180° sieve inserts can be replaced as desired in the corresponding retaining frames.

- ⇒ Remove the front of the corresponding retaining frame as described in Chapter "Cleaning the Grinding Set".
- Remove the existing sieve insert and replace it by the desired sieve insert instead.
- ⇒ Ensure the **correct orientation** when inserting sieve inserts with **trapezoidal perforation!**The direction of the arrow on the sieve insert must correspond to the direction of the arrow on the retaining frame! Both indicating the direction of rotation of the rotor (counterclockwise).
- ⇒ Screw the retaining frame back together.



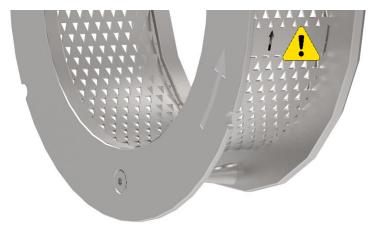


Fig. 22: Correct orientation of sieve inserts with trapezoidal perforation

# 6.8 Removing the Grinding Set

The removal of the grinding set is preferably carried out in the following sequence:

- 1. Rotor
- 2. Retaining frame
- 3. Grinding chamber cartridge

## 6.9 Mounting the Sample Receptacle

By the use of the textile filter hose (L), or a ring filter that is available as an optional accessory, attached between the discharge flange (AF) and the collecting receptacle (M), the airflow generated by the rotating rotor can be dissipated and the material throughput can be accelerated.

**NOTICE** If the collecting receptacle is installed without the filter hose or ring filter, it is to be expected that dust will be emitted out of the feed hopper (**F**). Therefore, never operate the SR 300 without the filter hose or ring filter!

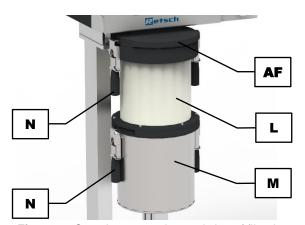
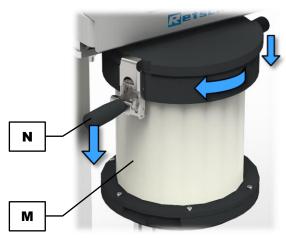


Fig. 23: Sample receptacle consisting of filter hose and collecting receptacle

- ⇒ Use the two horizontally unfolded clamping levers (**N**) to lift up the filter hose (**L**).
- ⇒ Position the filter hose (L) flush over the groove (AF1) of the discharge flange.
- ⇒ Turn the filter hose (L) clockwise until the clamping lever is situated over the clamping edge (AF2).
- $\Rightarrow$  Press the clamping levers (**N**) downwards with the open palm to clamp the filter hose (**L**).





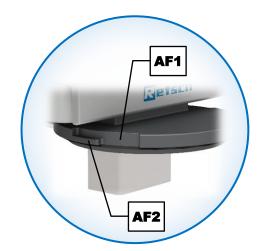


Fig. 24: Clamping the filter hose

- ⇒ Use the two horizontally unfolded clamping levers (N) to lift up the collecting receptacle (M).
- ⇒ Position the collecting receptacle (M) flush over the groove of the filter hose (L).
- ⇒ Turn the collecting receptacle (M) clockwise until the clamping lever is situated over the clamping edge.
- ⇒ Press the clamping levers (N) downwards with the open palm to clamp the collecting receptacle (M).

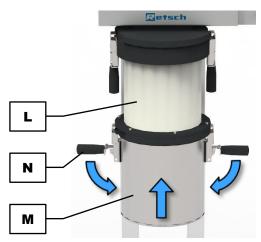


Fig. 25: Clamping the collecting receptacle

# **A** CAUTION

# **Danger of crushing**

Clamping the clamping levers

- If the clamping levers are completely enclosed by the hand, the fingers can be pinched.
- Do not clasp the clamping levers completely with the fingers during clamping.
- . Press down the clamping levers only with the palm.







### 7 Controlling the Device

### 7.1 Operating Controls, Display and Functions

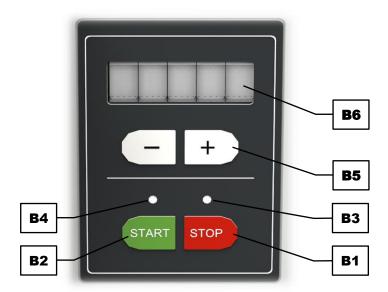


Fig. 26: Operating controls and functions

Element	Description	Function
B1	STOP button	Stops the grinding process
B2	START button	Starts the grinding process
В3	Status LED STOP	Lights up when STOP is pressed
В4	Status LED START	Lights up when START is pressed
В5	Rotation speed setting	Decreases or increases the rotation speed by pressing the "-" or "+" button, respectively in the range of 3 000 to 10 000 revolutions per minute
В6	Display	Displays the rotation speed as well as notifications

#### 7.2 Manual Mode

#### 7.2.1 Setting the Rotation Speed

The rotation speed can be set between 3 000 and 10 000 revolutions per minute (rpm) using the "+" and "-" buttons (**B5**).

- ⇒ Press the "+" button to increase the rotation speed in steps of 500 rpm.
- ⇒ Press the "–" button to decrease the rotation speed in steps of 500 rpm.

The rotation speed can also be changed during operation by pressing the "+" or "-" button (**B5**). An exceeding or falling below of 10 000 rpm or 3 000 rpm respectively is not possible.

#### 7.2.2 Start Process

⇒ Press the START button (B2) to start the grinding process.

The status LED (**B4**) lights up, the electrical locking mechanism is automatically activated and the rotor starts running.



**NOTICE** Always start the grinding process fist, before start feeding the sample material!

#### 7.2.3 Feeding Sample Material

**NOTICE** Start the SR 300 first, before start feeding the sample material!

The maximum feed grain size must not exceed 25 mm.

- ⇒ Fill the sample material slowly and continuously into the feed hopper (**F**) while the device is running.
- ⇒ Pay attention to the motor noise during the filling. If the rotor speed changes audibly due to excessive sample amount, immediately reduce the supply of the sample material.
- ⇒ During the grinding process, pay attention to the amount of grinding material in the collecting receptacle (**M**). The collecting receptacle must be emptied as soon as the filling level reaches 90 % of its volume.
- Tor the batch-wise or continuous operation, the SR 300 can be retrofitted with a 30 I collecting receptacle and the <u>vibratory feeder DR 100</u>, both available as optional <u>accessories</u>.

#### **NOTICE**

N14.0003

#### Damage to mechanical components

Rotor blocking

- Due to the high pulling power of the rotor, blockages can occur when feeding large-grained, harder sample material.
- Switch off the device immediately after blocking and remove the blocking grinding material.
- Reduce the feed of the sample material into the feed hopper.
- Fill larger and harder sample material slowly and gradually into the feed hopper.

#### 7.2.4 Stop Process

⇒ Press the stop button (B1) to stop the grinding process.

The status LED (**B3**) lights up, the rotor idles and the electric locking mechanism is automatically unlocked as soon as the rotor has come to a complete halt.

**NOTICE** Stop the grinding process only when there is no more sample material left in the grinding chamber!



### 8 Error Messages and Information Notes

### 8.1 Error Messages

Error messages inform the user about detected device or programme errors. In the event of an error message, a fault has occurred, in which the operation of the device or the programme is automatically interrupted. Such faults must be resolved before next startup.

Error code	Description	Measures
E10	Drive overload	⇒ Switch off the main switch and wait for 30 s
		before switching on again.
		⇒ If the error persists, contact service.
E11	Failure drive / motor	⇒ Switch off the main switch and wait for 30 s
		before switching on again.
		⇒ If the error persists, contact service.
E20	Failure main board	⇒ Switch off the main switch and wait for 30 s
		before switching on again.
		⇒ If the error persists, contact service.
E22	Failure keypad	⇒ Switch off the main switch and wait for 30 s
		before switching on again.
		⇒ If the error persists, contact service.
E26	Failure frequency	⇒ Switch off the main switch and wait for 30 s
	converter	before switching on again.
		⇒ If the error persists, contact service.
E41	Failure speed sensor	⇒ Switch off the main switch and wait for 30 s
		before switching on again.
		⇒ If the error persists, contact service.
E50	Failure safety circuit	⇒ Switch off the main switch and wait for 30 s
		before switching on again.
		⇒ If the error persists, contact service.
E51	Failure safety switch	⇒ Switch off the main switch and wait for 30 s
	(locking mechanism)	before switching on again.
		⇒ If the error persists, contact service.
E80	Failure interface	⇒ Switch off the main switch and wait for 30 s
		before switching on again.
		⇒ If the error persists, contact service.

#### 8.2 Information Notes

Notices inform the user on specific device or programme processes. The operation of the device or programme may be interrupted briefly, but there is no fault. The information notice must be acknowledged by the user to continue the process. Information notices provide additional information for the user as an aid, but do not represent any device or programme errors.

Notice code	Description	Measures
H10	Allow drive to cool down	⇒ Stop the grinding process.
		⇒ Allow the device to cool down.
H41	Close grinding chamber	⇒ Close the door.
H42	Open and close door	⇒ Open the door.



### 9 Return for Service and Maintenance



Fig. 27: Return form

The acceptance of devices and accessories of the Retsch GmbH for repair, maintenance or calibration can only be effected, if the return form including the decontamination declaration service has been correctly and fully completed.

- Download the return form located in the download section "Miscellaneous" on the Retsch GmbH homepage (<a href="http://www.retsch.com/downloads/miscellaneous/">http://www.retsch.com/downloads/miscellaneous/</a>).
- ⇒ When returning a device, attach the return form to the outside of the packaging.

In order to eliminate any health risk to the service technicians, Retsch GmbH reserves the right to refuse the acceptance and to return the respective delivery at the expense of the sender.



### 10 Cleaning, Wear and Maintenance

### A

### **CAUTION**

C14.0013

#### Personal injury

Improper repairs

- This manual does not contain any repair instructions.
- For safety reasons, repairs may only be carried out by Retsch GmbH or an authorised representative or by qualified service technicians.

#### 10.1 Cleaning



#### **WARNING**

W4.0003

#### Danger to life through electric shock

Cleaning with water

- An electric shock can cause burns, cardiac arrhythmia, respiratory arrest, as well as cardiac arrest.
- The power cable must be unplugged before cleaning the device.
- Use a cloth dampened with water for cleaning.
- Do not clean the device under running water!





### A

#### **CAUTION**

C15.0031

#### Risk of injury

Cleaning with compressed air

- When using compressed air for cleaning purposes dust and remnant of the sample material can be flung around and injure eyes.
- · Always wear safety glasses when cleaning with compressed air.
- . Observe the material safety data sheets of the sample material.





#### **NOTICE**

N15.0009

#### Damage to the housing and device

Use of organic solvents

- Organic solvents may damage plastic parts and the coating.
- The use of organic solvents is not permitted.
- Clean the housing of the device with a damp cloth and if necessary, with a household cleaning agent. Pay attention that no water or cleaning agent enters the interior of the device.
- ⇒ Clean the grinding chamber and the motor shaft (**MW**) with a brush or paint brush and vacuum the loosened material residues with a vacuum cleaner.
- ⇒ Alternatively, the grinding chamber can also be cleaned with compressed air.



### 10.1.1 Cleaning the Grinding Set

The cleaning of the grinding set consisting of the grinding chamber cartridge (**MK**), the retaining frame with the sieve insert (**SR**) and the rotor (**RO**) should be carried out regularly.

⇒ Remove the grinding set from the grinding chamber (→ Chapter "Removing the Grinding Set")

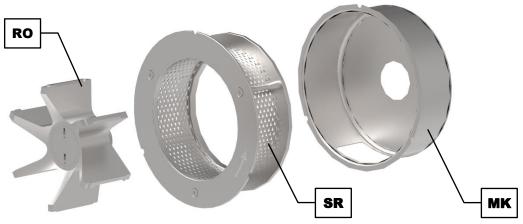


Fig. 28: Grinding set

#### Cleaning the grinding chamber cartridge:

The grinding chamber cartridge (**MK**) can be cleaned with alcohol, petrol or normal household detergent. The grinding chamber cartridge is also dishwasher suitable.

#### Cleaning the retaining frame:

For thorough cleaning of the retaining frame for 360° sieve inserts, it is recommended to separate the sieve insert (**SB**) from the retaining frame (**SR**).

- ⇒ Remove the three M5 hexagon socket countersunk head screws.
- ⇒ Remove the front retaining frame to remove the 360° sieve insert.

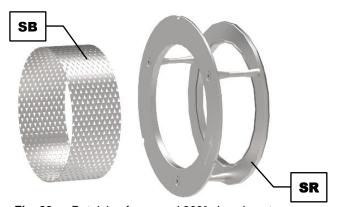


Fig. 29: Retaining frame and 360° sieve insert



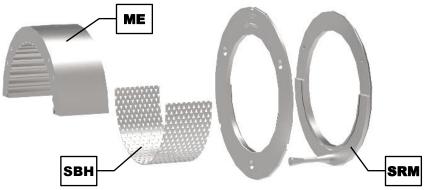


Fig. 30: Retaining frame, grinding insert and 180° sieve insert

For thorough cleaning of the retaining frame for 180° sieve inserts, it is recommended to separate the sieve insert (**SBH**), as well as the grinding insert (**ME**) from the retaining frame (**SRM**).

- ⇒ Remove the three M5 hexagon socket countersunk head screws.
- ⇒ Remove the front retaining frame to remove the 180° sieve insert.
- Additionally remove the two M5 hexagon socket countersunk head screws that connect the grinding insert and the rear retaining frame.

The retaining frames (**SR**), (**SRM**), the sieve inserts (**SB**), (**SBH**), as well as the grinding insert (**ME**) can be cleaned with alcohol, petrol or normal household detergent. Sieve inserts (**SB**), (**SBH**) can be easily and effectively cleaned dry or wet with a hand brush with plastic bristles. All parts are also dishwasher suitable.

Sieve inserts (**SB**), (**SBH**) with finer mesh sizes can also be cleaned in an ultrasonic cleaning-bath. As cleaning agent, water together with a standard surfactant is recommended. The cleaning in the ultrasonic bath usually takes two to three minutes. After that the sieve inserts are thoroughly rinsed with water and dried. The cleaning with strong bases or acids is generally not recommended.

Drying ovens of various sizes can be used for drying sieve inserts.

Additional information concerning ultrasonic cleaning-baths and drying ovens can be found on the Retsch GmbH homepage (http://www.retsch.com).

#### Cleaning the rotor:

The rotor (**RO**) can be cleaned with alcohol, petrol or normal household detergent. The rotor is also dishwasher suitable.

Make sure that the hub (RO1) is thoroughly cleaned and no material residues are left inside.



### 10.1.2 Cleaning the Feed Hopper

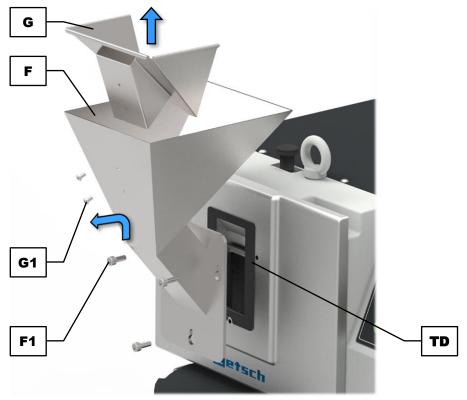


Fig. 31: Cleaning the feed hopper

- ⇒ Unscrew the two M4 oval-head screws (G1).
- ⇒ Remove the splashback protection (**G**).
- ⇒ Loosen the three M6 hexagon socket head screws (F1).
- ⇒ Pull the feed hopper (**F**) slightly upwards, then remove it to the front.
- ➡ Clean the component parts with alcohol, petrol or normal household detergent. The feed hopper (F) is also dishwasher suitable.
- ⇒ Check the sealing gasket (**TD**) on the door for its intactness and cleanliness before reinstalling the feed hopper (**F**) and the splashback protection (**G**) after cleaning.

## **WARNING**

W5.0011

### Serious personal injury

Intervention in rotating rotor



- Unintentional intervention in the grinding chamber and the rotating rotor can cause serious hand injuries.
- Never operate the device without mounted splashback protection, or, if removable, without mounted feed hopper.

#### 10.2 Wear

The grinding tools may become worn, depending on the frequency of the grinding operation and the sample material. The rotor and the grinding sets should be regularly checked for wear and replaced if necessary.



Likewise, all existing sealing gaskets (of grinding tools and in the device) should be checked for wear on a regularly basis and replaced if necessary.

#### 10.3 Maintenance

The SR 300 is largely maintenance-free.

With each cleaning, it is recommended to additionally lubricate the feather key (**PF**) on the motor shaft (**MW**) with a few drops of machine oil.

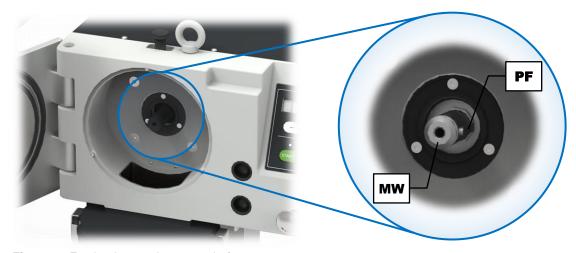


Fig. 32: Feather key on the motor shaft

In order to ensure the operational safety of the device, the rollers (K1) of the locking mechanism (K) should be checked monthly for free movement and, if necessary, re-oiled with machine oil. At the same time, the magnets (K2) should be thoroughly cleaned.

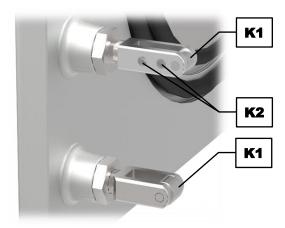


Fig. 33: Locking mechanism



#### 11 Accessories

Information on available accessories as well as the respective manuals are accessible directly on the Retsch GmbH homepage (http://www.retsch.com) under the heading "Downloads" of the device.

Information on wear parts and small accessories can be found in the Retsch GmbH general catalogue also available on the homepage.

In case of any questions concerning spare parts please contact the Retsch GmbH representative in your country, or Retsch GmbH directly.

### 11.1 Vibratory Feeder DR 100



Fig. 34: SR 300 with vibratory feeder DR 100

When feeding larger quantities, it is generally recommended to use the vibratory feeder DR 100 for uniform conveyance of the feed material. This mainly avoids an unnecessary load on the grinding set and ensures reproducibly exact results, as well as the economical use of the downstream device.

**NOTICE** Observe the supplied manual before installing and operating the vibratory feeder DR 100.



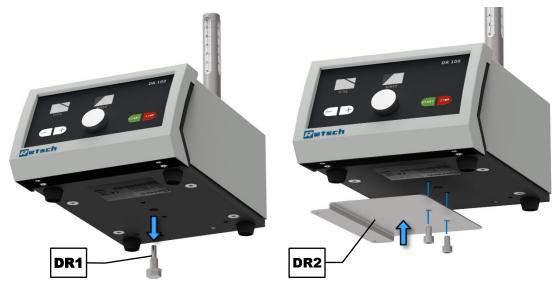


Fig. 35: Removing the transportation lock (left) and mounting the angle plate (right)

- ⇒ Remove the transportation lock (**DR1**).
- ⇒ Mount the angle plate (**DR2**) to the bottom of the DR 100 with the two M6 hexagon socket head screws supplied, so that it protrudes under the vibratory feeder on the left hand side.

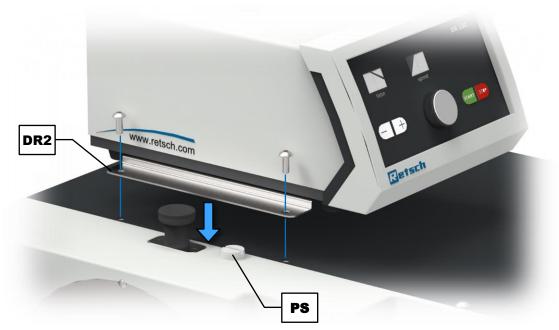


Fig. 36: Mounting the vibratory feeder DR 100

- ⇒ If not already done, replace the transportation lug (**A**) on the SR 300 with the supplied plastic screw (**PS**).
- ⇒ Loosen the two M5 oval-head screws at the housing of the SR 300 and use them to screw the angle plate (**DR2**) to the SR 300.
- ⇒ Mount the feed chute holder, the feed chute, the hopper holder and the hopper as described in the separate manual of the vibratory feeder DR 100.





Fig. 37: Rear side of the vibratory feeder DR 100

- ⇒ Plug the power cable into the mains connection on the rear.
- ⇒ Set the operating mode selector switch (**DR3**) to "Standard". A control via "Extern" (external) is not possible in conjunction with the SR 300.
- ⇒ Turn on the vibratory feeder DR 100 via the rear mains switch (**DR4**).
- ⇒ For detailed information on the controls, please refer to the manual of the vibratory feeder DR 100.

### 11.2 Cyclone Separator

## **WARNING**

#### Serious personal injury

Intervention in rotating rotor

- Unintentional intervention in the grinding chamber and the rotating rotor can sever fingers and cause serious hand injuries.
- The power cable must be unplugged before removing the discharge flange.
- Never operate the device without mounted discharge flange.







**Fig. 38:** SR 300 and cyclone separator with 5 I collecting receptacle (left) and 30 I collecting receptacle (right)

When grinding light sample material, the SR 300 can be operated with a cyclone separator, so that even light feed material or small quantities can be easily processed.

Depending on the amount of sample to be processed, the cyclone separator can be fitted with a 5 I or 30 I collecting receptacle.

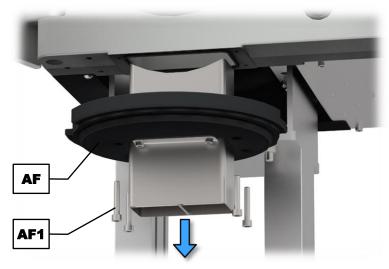


Fig. 39: Removing the discharge flange

- ⇒ Disconnect the device from the mains.
- ⇒ Loosen the four M6x35 hexagon socket head screws (**AF1**).
- ⇒ Remove the discharge flange (**AF**).



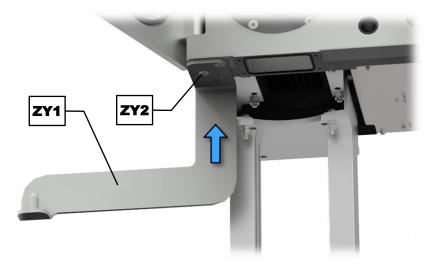


Fig. 40: Installing the mounting bracket

- ⇒ Use the three supplied M8 hexagon socket countersunk head screws (**ZY2**) to bolt the mounting bracket (**ZY1**) of the cyclone separator to the housing of the SR 300.
- ① Once installed, the mounting bracket (**ZY1**) does not need to be removed again for the operation without cyclone separator. The discharge flange (**AF**) can also be reassembled safely with the mounting bracket installed.


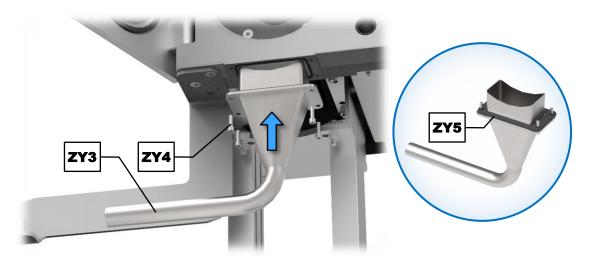


Fig. 41: Installing the cyclone adapter

- ⇒ Fasten the cyclone adapter (**ZY3**) to the housing of the SR 300 with the four supplied M6x20 hexagon socket head screws (**ZY4**).
- ⇒ Make sure that the sealing gasket (**ZY5**) rests correctly on the cyclone adapter.



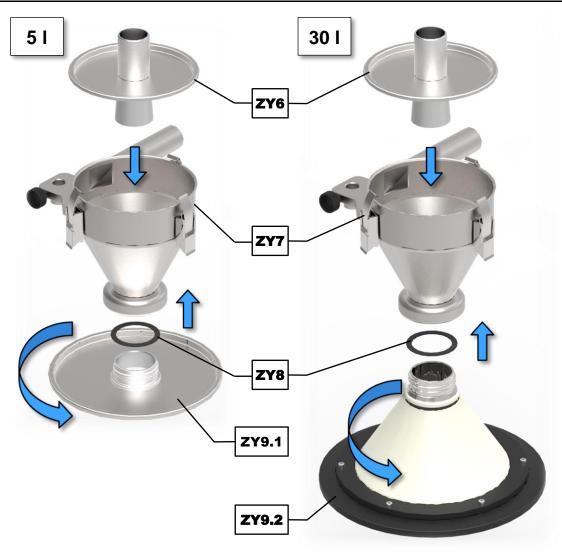

**Fig. 42:** Assembly of the cyclone funnel for the 5 I collecting receptacle (left) and the 30 I collecting receptacle (right)

- ⇒ Depending on the type of the cyclone separator, either screw the cover (**ZY9.1**) of the 5 I collecting receptacle together with the sealing gasket (**ZY8**) into the socket of the cyclone funnel (**ZY7**), or, in case of the 30 I collecting receptacle, screw the filter hose (**ZY9.2**) together with the sealing gasket (**ZY8**) into the socket of the cyclone funnel (**ZY7**).
- ⇒ Open the three snap locks on the cyclone funnel (**ZY7**), place the funnel cover (**ZY6**) onto the cyclone funnel and close the three snap locks again.



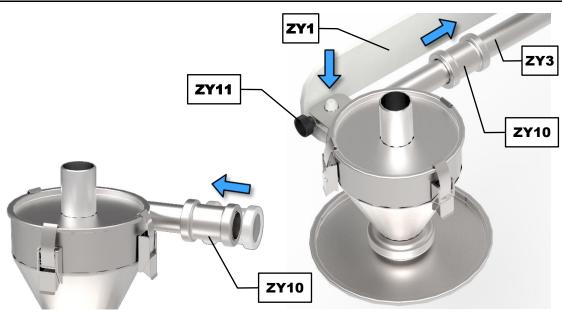


Fig. 43: Placing the coupling (left) and mounting the cyclone funnel (right)

- ⇒ Slide the coupling (**ZY10**) completely over the tube of the cyclone funnel.
- ⇒ Mount the assembled cyclone funnel into the mounting bracket (**ZY1**) and align it to the cyclone adapter (**ZY3**).
- ⇒ Slide the coupling (**ZY10**) over the cyclone adapter (**ZY3**) until it is situated half on the tube of the cyclone funnel and half on the cyclone adapter.
- ⇒ Secure the assembled cyclone funnel by means of the knurled head screw (ZY11).

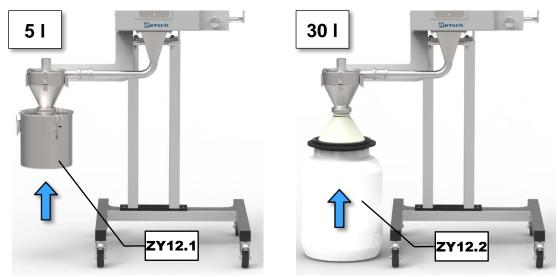


Fig. 44: Inserting the 5 I (left) and 30 I collecting receptacle (right)

⇒ Depending on the type of the cyclone separator, clamp the 5 I collecting receptacle (**ZY12.1**) to the cover (**ZY9.1**) by means of the three snap locks, or place the 30 I collecting receptacle (**ZY12.2**) underneath the filter hose (**ZY9.2**) so that its flange rests on the opening of the collecting receptacle.





Fig. 45: Connecting the industrial vacuum cleaner

- ⇒ Insert the connection of the industrial vacuum cleaner (**ZY13**) into the upper opening of the cyclone funnel.
- ⇒ Always switch on the industrial vacuum cleaner first, before starting the grinding process.

**A** CAUTION Read the manual of the industrial vacuum cleaner before commissioning.



### 12 Disposal

In the case of a disposal, the respective statutory requirements must be observed. In the following, information on the disposal of electrical and electronic devices in the European Community are given.

Within the European Community the disposal of electrically operated devices is regulated by national provisions that are based on the EU Directive 2012/19/EU on Waste Electrical and Electronic Equipment (WEEE).

Accordingly, all devices supplied after August 13<sup>th</sup> 2005 in the business-to-business area, to which this product is classified, may no longer be disposed of with municipal or household waste. To document this, the devices are provided with the disposal label.

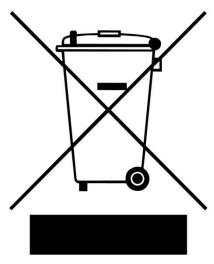


Fig. 46: Disposal label

Since the disposal regulations worldwide and also within the EU may differ from country to country, the supplier of the device should be consulted directly in case of need.

This labelling obligation is applied in Germany since March 23<sup>rd</sup> 2006. From this date on, the manufacturer must provide an adequate possibility of returning all devices delivered since August 13<sup>th</sup> 2005. For all devices delivered before August 13<sup>th</sup> 2005 the end user is responsible for the proper disposal.



# 13 Index

1	E
180° sieve insert13, 43	Elec
3	Elec
360° sieve insert13, 42	EMC
A	Eme
••	Emi
Accessory 46	Equi Erro
base frame20	E <sup>2</sup>
Action instructions7	E,
Ambient temperature15, 16	E
Amperage17	E
Application-technical information6	E2 E4
В	E:
Back28	E:
Back view28	E
Bar code17	Erro
Base frame20	Expl
C	Exte
Calibration40	F
Capacity17	Feat
Catch mechanism28	Fee
CE marking	Fee
Clamping lever	cle
Cleaning	Fee
Collecting receptacle28, 35	Fee
Complaints14	Filte Fina
Condensation15	First
Confirmation form for the managing operator10	Fred
Convigation 6	Fror
Copyright6 Cutting injuries32	Fron
Cyclone separator	Fund
D	Fuse Fuse
Degree of protection11	
Depth12	G
base13	Gen
with open door13	Gen Grin
Device	cle
close29	in
open29	Grin
Device designation17	clo
Dimensions12	Grin
Disclaimer	Grin
Disconnection from the mains	Grin
Display	cle
label54	in
regulations54	re
Disposal label17	Н
Distance rotor12	Hea

Electrical connection	
Electromagnetic compatibility EMC	
Emergency unlocking	
Emissions	
Equivalent continuous sound level 11,	12
Error E10	30
E11	
E20	
E22	
E26	
E41 E50	
E51	
E80	
Error messages	39
Explanations of the safety instructions	7
External fuse	16
F	
Feather key	
Feed grain size	
Feed hopper	
cleaning	
Feed volume Feeding sample material	
Filter hose28,	35
Final fineness	
First commissioning	
Frequency	
Front view	
Functions	
Fuse strength	
Fuse type	17
G	
General catalogue	46
General safety instructions	8
Grinding chamber cartridge	40
cleaninginsert	
Grinding insert	
cleaning	
Grinding materials	
Grinding noises	
Grinding set	42
cleaning	
insertremove	
	ა၁
H	
Hearing damage	
Height	12



Humidity16	Repair	9, 40, 41
ı	Repair instructions	
Information mate	Required floor space	13
Information note	Retaining frame	
H1039	cleaning	
H4130, 39 H4230, 39	for 180° sieve insert	
•	for 360° sieve insert	
Information notes39	insert	34
Installation14	Return	14
Installation height15	for service and maintenance	40
Installation site	Return device	54
conditions15	Return form	
L	Revision status	
L <sub>eq</sub> 11, 12	Ring filter	
Lifting the device	Rotation speed	
by persons15	setting	37
with hoist15	Rotor	
Location requirements13	cleaning	
Locking lever28, 29	insert	
Locking mechanism11, 28, 45		
electrical	Round perforation	13
mechanical	S	
•	Safety instruction	7
Locking pin28, 29	caution	
Long-term operation26	danger	
М	notice	
Mains connection29	warning	
Mains frequency17	Safety manager	
Mains supply16	Sample receptacle	
Mains switch29	mount	35
Maintenance 10, 40, 41, 45		
Manual6, 8, 10	Serial number	
Manual mode37	Service address	9
Manufacturer's address17	Setting up	4.0
Motor rotation speed12	as table-top device	
N '	on base frame	
	Sieve diameter	
Notes on the Manual6	Sieve insert	
0	change	
Opening30	cleaning	
Operating controls28, 37	cleaning of fine mesh sizes	
Operating instructions10	drying	
Operating the device25	Signs	
P	Small accessories	
r	Sound level	
Packaging14	Sound parameters	
Part number17	Spare parts	
Power version17	Splashback protection	
Principle of operation26	START button	
Process	Status LED	
start37	STOP button	
stop38	Switching on / off	
Protective equipment11	Symbols	7
R	Т	
	Target group	g
Range of application of the device26	Technical data	
Rated power	Temperature fluctuations	
Receptacle volume	Temperature range	
Relative humidity	Temporary storage	





Transport Transport shakles remove	
Transportation damage	.29 .13
Type plate16, 17, description	.17
UKCA marking Use of the device for the intended purpose	
VibrationsVibratory feeder DR 100	

Views of the instrument27	7
Voltage16	ô
W	
Warranty claims 9, 14	4
Wear 32, 41, 44	4
Wear parts 46	ဝ
Weight 12, 15, 17	
Width12	2
base13	
with open door13	3
Working space13	3
Workplace related emission level 11, 12	
Υ	
Year of production17	7



### **EU Declaration of Conformity**

Translation

### **ROTOR BEATER MILL**

SR 300 | 20.752.xxxx

#### **EU DECLARATION OF CONFORMITY**

We, represented by the undersigned, hereby declare that the above device complies with the following directives and harmonised standards:

#### Machinery directive 2006/42/EC

Applied standards, in particular:

DIN EN ISO 12100:2011 Machine Safety - General Design Principles

DIN EN 13683:2013 Garden equipment – Integrally powered shredders/chippers – Safety
DIN EN 61010-1 Safety Regulations for Electrical Measurement, Control, Regulation and

Laboratory Devices

#### Electromagnetic compatibility 2014/30/EU

Applied standards, in particular:

EN 55011:2009 Industrial, scientific and medical equipment - Radio-frequency disturbance

+A1:2010 characteristics - Limits and methods of measurement

DIN EN 61326-1:2013 Electrical equipment for measurement, control and laboratory use - EMC

requirements

#### Restriction of hazardous substances (RoHS) 2011/65/EU

Applied standards, in particular:

DIN EN IEC 63000:2019 Technical documentation for the assessment of electrical and electronic

products with respect to the restriction of hazardous substances

#### Authorised person for compilation of the technical documentation:

Julia Kürten (Technical Documentation)

Furthermore, we declare that the relevant technical documentation for the above device has been prepared in accordance with Annex VII Part A of the Machinery Directive and we undertake to submit this documentation to the market surveillance authorities on request.

In the event of a modification of the device not agreed on by Retsch GmbH, as well as the use of non-approved spare parts or accessories, this declaration loses its validity.

Retsch GmbH Haan, 01/2023

Dr. Kevin Schmitz, Head of Development

part of VERDER

© Retsch GmbH · Retsch-Allee 1-5 · 42781 Haan · Germany www.retsch.com · E-mail: info@retsch.com · Phone: +49 2104 2333-100





# Copyright

© Copyright by Retsch GmbH Retsch-Allee 1-5 42781 Haan Germany