

EasyMax 402

Personal Synthesis Workstation



METTLER TOLEDO

Table of Contents

1	Introduction	3
1.1	Scope of delivery	3
1.2	Check on arrival	4
2	Safety Information	5
2.1	Definition of signal warnings and symbols.....	5
2.2	Intended use	5
2.3	Product-specific safety.....	5
3	Overview	8
4	Installation	9
4.1	Installation requirements	9
4.2	Unpacking and transporting device	9
4.3	Connecting power to the device	9
4.4	Connecting a flowmeter	10
4.5	Installing instrument cooling	10
4.5.1	Connecting in-house coolant supply.....	11
4.5.2	Connecting cryostat cooling	11
4.6	Connecting instrument purge.....	12
4.7	Connecting touchscreen to EasyMax	12
4.8	Assembling the stirrer	12
4.9	Installing a Tr sensor.....	13
4.10	Installing a two-piece 400 mL reactor	13
4.11	Turn on Device	14
5	Operation	15
5.1	Select Reactor Type	15
5.2	Change Safety Settings	16
5.2.1	Change Safety Temperature (T safe)	16
5.2.2	Change Reaction Temperature limits (Tr)	17
5.2.3	Change Range of Jacket Temperature (Tj)	17
5.2.4	Change T diff max	17
5.2.5	Change Rsafe	18
5.2.6	Change Rmax.....	18
5.3	Start an Experiment	19
5.4	Change Stirrer Speed	19
5.5	Change Tj	19
5.6	Change Tr	20
5.7	End an Experiment.....	20
6	Maintenance	22
6.1	Update Firmware.....	22
6.2	Checking the Reactor.....	22
6.3	Cleaning the Instrument	22
6.4	Disposal	22
7	Technical Data	23
7.1	Thermostat	24
7.2	Reactors	25
7.3	Stirrer	25
7.4	Cooling	26
7.5	Purge gas	26

1 Introduction

The METTLER TOLEDO EasyMax 402 is an easy-to-use personal synthesis workstation that allows you to run two reactions simultaneously in an accurate and reproducible way. The EasyMax 402 has two reactor positions that can be controlled independently designed for a variety of reactors and vials. The unique cooling principle allows to run experiments from -40 °C to 180 °C without the use of oil or ice baths, or a bulky cryostat. Additionally, control your experiments with an intuitive touchscreen.








The main features of the EasyMax 402 are:

- Two independently controlled reactor zones (for 100 mL and/or 400 mL one or two-piece reactors with glass or PTFE covers)
- Overhead stirring
- No oil or ice baths
- Easy touchscreen control
- Full compatibility with selected METTLER TOLEDO devices, including the DU SP-50, ECB, EasySampler and the SevenExcellence pH meter

Please also read the Operating Instructions for the full scope of functionalities of the device. The Operating Instructions can be found on the USB stick.

1.1 Scope of delivery

The following items are included in the EasyMax 402 Basic and EasyMax 402 Advanced thermostat set:

Order number		Description	Quantity
		EasyMax 402 thermostat	1
51161883		TFT touchscreen 7", 1 m cable	1
11132570		Protective cover for touchscreen	1
51191125		PVC hose, soft, for reflux condenser, 5 m	1
51161187		PVC industrial hose for coolant, 15 bar, 2.5 m	2
51192239		PVC industrial hose for purge gas, 18 bar, 2 m	1
51161186		PVC hose for purge gas, 2 m	2
51191373		Y-piece for gas tubing	2

51191916		Reducing connector for purge gas tubing	1
51161827		Flowmeter set	1
51190324		Quick connect coupling for purge gas inlet	4
51192126		Hose clamp for PVC tube	8
51191915		Flow indicator for coolant	1
51191914		Knurled screw, M6 x 10	3
51191235		Identification clips for adhesive labels	10
		Set of adhesive labels	1
30356209		USB Stick RXE/CSS with documents	1
		User Manual	1

Additional items shipped with EasyMax 402 Advanced

	iControl software CD package	1
	iC Data Center voucher for EasyMax/OptiMax (voucher only), one year	1
	iControl Office License (license sheet only)	1

If an item is missing, please contact your local support team.

1.2 Check on arrival

Check the following conditions once the package has arrived:

- The package is in good condition.
- The contents show no signs of damage (e.g. broken covers, scratches, etc.)
- The content is complete (see [Scope of delivery ► Page 3]).

If any one of these condition is not fulfilled, please contact your local support team.

2 Safety Information

This thermostat has been tested for the intended purposes described in this document. However, this does not absolve you from the responsibility of performing your own tests of the product supplied by us regarding its suitability for the methods and purposes you intend to use it for. You should therefore observe the following safety measures.







We, Mettler-Toledo GmbH, accept no liability whatsoever if you do not observe the following rules and safety notes for safe operation of the thermostat.

2.1 Definition of signal warnings and symbols

Safety notes are indicated by signal words and warning symbols and contain warnings and information about safety issues. Ignoring safety notes can lead to personal injury, damage to the instrument, malfunctions and erroneous results.

WARNING	A hazardous situation with medium risk, possibly resulting in death or severe injury if not avoided.
CAUTION	A hazardous situation with low risk, resulting in minor or moderate injury if not avoided.
NOTICE	A hazardous situation with low risk, resulting in damage to the instrument, other material damage, malfunctions and erroneous results, or loss of data.
Note	(no symbol) for useful information about the product.

Meaning of safety symbols

	Electrical Hazard		Explosion		Burns / Hot Surface
	Rotating parts		Heavy load		General note

2.2 Intended use

The METTLER TOLEDO EasyMax 402 is a reactor system for performing parallel synthesis, with 100-mL and 400-mL glass reactors.

The device is designed to be used in a laboratory environment and operated in a fume hood. All users should be trained to work in a laboratory and with this device.

Always operate and use your device in accordance with the instructions contained in this manual; use it only together with equipment specified in this documentation.

Any other type of use and operation beyond the limits of these technical specifications without the written consent from Mettler-Toledo GmbH is considered as not intended.

2.3 Product-specific safety



WARNING

Risk of electric shock

- 1 Make sure to plug the power cable supplied into a power supply outlet that is grounded. A technical fault could otherwise result in serious injury or death.
- 2 Only use the METTLER TOLEDO power supply cable and AC power adapter designed for your instrument.



WARNING

Power failure

A power failure can lead to explosion with possibly fatal consequences.

- Implement appropriate measures like an uninterruptible power supply (UPS).



WARNING

Risk of explosion with critical reactions

Performing critical reactions could lead to explosions.

- Perform a safety analysis before starting an experiment with high hazardous potential for example by using a Differential Scanning Calorimeter.



WARNING

Risk of explosion due to damaged reactors

Explosion of a reactor could cause serious injury.

- Check the reactor before each use for damage (scratches, formation of cracks).



CAUTION

Hot parts when working above 50 °C

Touching hot parts can cause burns.

- Do not touch the cover plate of the device, the fixing ring, the reactor covers, attachments of the reactor or the overhead stirrer if you work above 50 °C.



CAUTION

Rotating parts of stirrer

Rotating parts of a running stirrer may lead to injuries.

- 1 Do not touch rotating parts of a stirrer.
- 2 Do not wear loose clothing and make sure jewellery and long hair do not get entangled in the stirrer.



NOTICE

Wrong coolant used

High chloride concentration or some additives in the coolant can lead to corrosion of the thermostat.

- 1 Do not use solutions of NaCl, CaCl₂ or DW-Therm.
- 2 Check compatibility with the wetted parts of the coolant system.



NOTICE

Wrong connection or disconnection of cables

A wrong connection or disconnection of cables during operation could lead to instrument damage.

- 1 Before switching the device on, connect the cables of stirrers and sensors to their respective inputs and outputs.
- 2 Do not disconnect the cables while the instrument is operating.



NOTICE

Condensation of atmospheric moisture

The condensation of atmospheric moisture can cause corrosion of the instrument.

- To prevent this, purge the instrument with dry air, nitrogen or argon.



NOTICE

Thermal shock

Glass parts of the instrument or the reactor could get damaged.

- Do not pour cold liquids into hot glassware and vice versa.



NOTICE

Risk of reactor breakage when sealed hermetically

Hermetical sealing could lead to pressure build-up when using gas or when the reactor is heated.

- Make sure venting is always possible.



NOTICE


Reactor breaking due to freezing

The reactor can break or get stuck in the reactor zone of the thermostat when atmospheric moisture or any liquids freeze on the outside of the reactor or in the reactor zone of the thermostat.

- 1 Make sure the reactor zone of the thermostat and the reactor itself are clean and dry before inserting.
- 2 Make sure that when cooling below 0 °C, there is enough purge to minimize condensation.


3 Overview

Front view

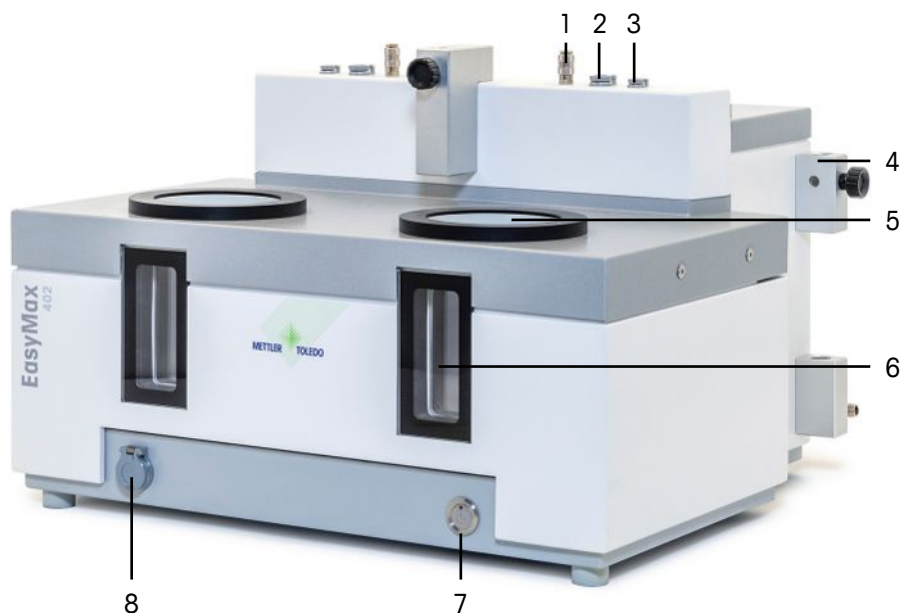

1	Reactor purge gas outlet (quick connect coupling)	2	Stirrer connection
3	Tr sensor port	4	Holder for lab bars
5	Opening for reactors	6	Window (with backlight)
7	Power button	8	USB Socket

Back view

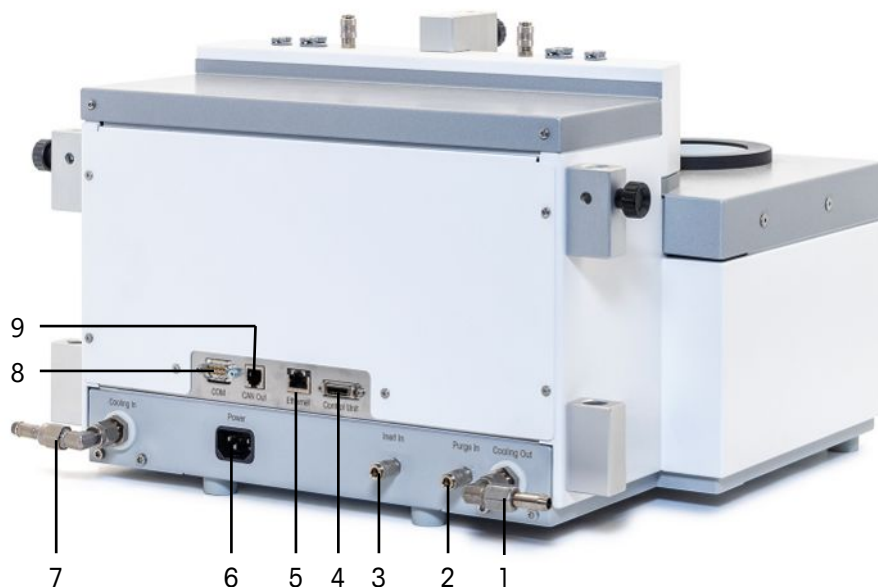

1	Cooling Out	2	Purge In connection
3	Inert In connection	4	Touchscreen connection
5	Ethernet connection	6	Power supply
7	Cooling In	8	RS232
9	CAN out		

4 Installation

4.1 Installation requirements

- The device should be installed in a fume hood.
- Make sure you install the device in accordance with the technical data.

Site requirements

The instrument has been developed for indoor operation in a well-ventilated area. Avoid the following environmental influences:

- Conditions outside of the ambient conditions specified in the technical data
- Powerful vibrations
- Direct sunlight
- Corrosive gas atmosphere
- Explosive atmosphere of gases, steam, fog, dust and flammable dust
- Powerful electric or magnetic fields

4.2 Unpacking and transporting device



CAUTION

Risk of injury due to heavy load

You can injure yourself by carrying the instrument alone.

- Never try to carry the instrument alone. At least two people are needed to carry the instrument.

Unpack the device

- 1 Grip the device under the base plate.
- 2 Lift the device up and out of the foam packing material.
- 3 Place the device on the lab bench.

Transport the device

- 1 Unplug the power adapter.
- 2 Disconnect the device properly from the cooling media.
- 3 Grip the device under the base plate.

4.3 Connecting power to the device



WARNING

Risk of electric shock

- 1 Make sure to plug the power cable supplied into a power supply outlet that is grounded. A technical fault could otherwise result in serious injury or death.
- 2 Only use the METTLER TOLEDO power supply cable and AC power adapter designed for your instrument.

- 1 The power supply connection is on the back side of the device.
- 2 Connect the instrument to the power supply using the included country-specific cable.
- 3 Insert the plug of the power cable in a grounded power outlet that is easily accessible.



4.4 Connecting a flowmeter




NOTICE

Damage of Flowmeter

The flowmeter is not resistant to any liquids other than water! Do not use it with other coolants.

- 1 Screw the flowmeter with the knurled screw onto the metal holder.
- 2 Insert the holder into the space between the lab bar holders.
- 3 Slide the lab bar through the lab bar holes and the holes of the metal holder.
- 4 Tighten the screw on the lab bar holder to fasten the lab bar.

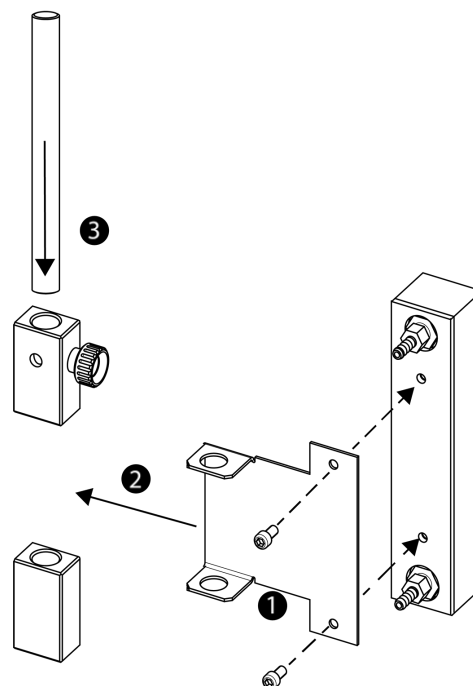

4.5 Installing instrument cooling

There are two possibilities to cool the instrument.



NOTICE

Damage of Flowmeter

The flowmeter is not resistant to any liquids other than water! Do not use it with other coolants.



NOTICE


Wrong coolant used

High chloride concentration or some additives in the coolant can lead to corrosion of the thermostat.

- 1 Do not use solutions of NaCl, CaCl₂ or DW-Therm.
- 2 Check compatibility with the wetted parts of the coolant system.

4.5.1 Connecting in-house coolant supply

The instrument needs a constant flow of the cooling medium.

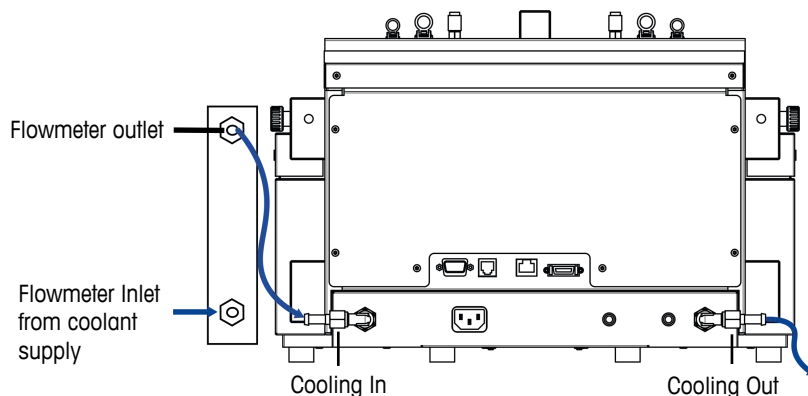

The instrument is delivered with two PVC industrial tubes. If you use the flowmeter, you need a third piece which can be cut away from one of the delivered tubes.

- 1 Push one piece of the PVC industrial hose (51161187) over the cooling inlet of the coolant flowmeter.
- 2 Secure it with a hose clamp.
- 3 Connect it to the coolant supply.
- 4 Push the cut piece of the PVC tube over the outlet of the coolant flowmeter.
- 5 Connect the other end to the elbow coupling of the **Cooling In** on the back side of the instrument.
- 6 Secure both connections with a hose clamp.
- 7 Push the other PVC tube over the elbow coupling of the **Cooling Out** on the back side of the instrument.
- 8 Secure it with a hose clamp.
- 9 Connect it to the cooling outlet of the fume hood.

4.5.2 Connecting cryostat cooling



NOTICE

Damage of Flowmeter


The flowmeter is not resistant to any liquids other than water! Do not use it with other coolants.

The instrument needs a constant flow of the cooling medium.

- 1 Screw the insulated connection tube over the **Cooling In** on the back of the instrument.
- 2 Use a wrench to tighten the connection to the instrument.
- 3 Screw the second insulated connection tube over the **Cooling Out** on the back of the instrument.
- 4 Use a wrench to tighten the connection to the instrument.

4.6 Connecting instrument purge

To prevent corrosion by condensed atmospheric moisture, purge the instrument with a dry gas, e.g. dry air, dry nitrogen or dry argon.

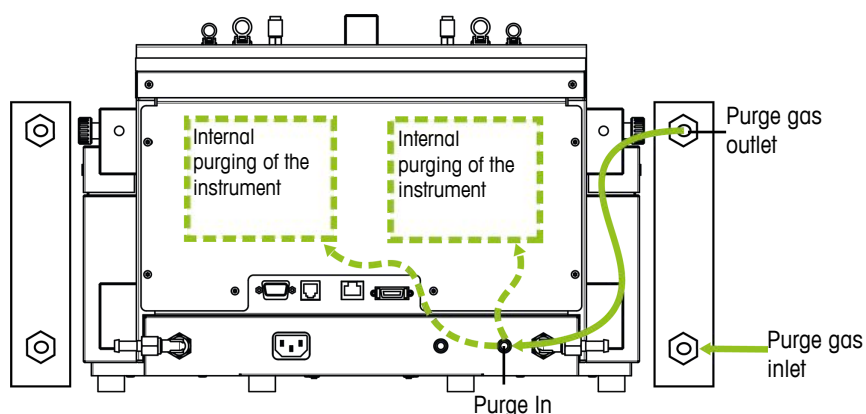

- 1 Install the quick connect coupling (51190324) with the PVC tubing (51161186) on the purge gas inlet connector.
- 2 Secure it with a hose clamp.
- 3 Connect the PVC tubing to the gas supply.

4.7 Connecting touchscreen to EasyMax

- 1 The touchscreen connection is on the back side of the instrument (see picture).
- 2 Connect the touchscreen cable to the socket.
- 3 Position the touchscreen so that it is always readable.



4.8 Assembling the stirrer

The overhead stirrer consists of the stirrer motor and a stirrer shaft that can be equipped with several stirrer elements (see Synthesis Workstation Catalog).



NOTICE


Glass breakage due to expanding material

The stirrer adapter is made of PEEK. This material can swell in combination with certain solvents, which may cause the glass to break.


- Use a PTFE stirrer adapter when working with these solvents.

- 1 Screw the pitched-blade element (7) onto the stirrer shaft (6).

- 2 Push the stirrer shaft (6) or the glass stirrer from below through the opening of the reactor cover.
- 3 Push the adapter (5) with the height adjustment unit (4) over the stirrer shaft and insert it into the tapered opening of the cover.
- 4 Place the PA washer (3) over the stirrer shaft onto the adjustment unit (4).
- 5 Push the lower coupling (2) onto the stirrer shaft.
- 6 Push the stirrer motor (1) over the adjustment unit (4) in such a way that it snaps into its three pins.
- 7 Check that the stirrer blade does not touch any inserts.
- 8 Secure the adapter on the reactor with a Keck clamp.

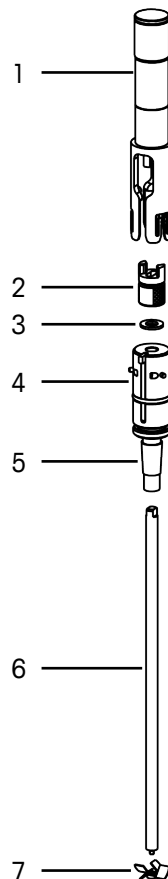

4.9 Installing a Tr sensor

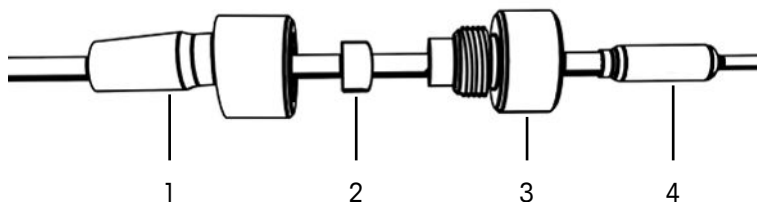

- 1 Slide a UNF 1/4" G 28 nut (2) over the sensor (4).
- 2 Slide a UNF 1/4" G 28 ferrule (3) over the sensor with the narrow end facing the nut.
- 3 Screw the nut lightly into the adapter (1) to press the ferrule into the nut.
- 4 Check that the Tr sensors does not touch the stirrer blades or other inserts.

Note Make sure the temperature sensor shows a sufficient immersion depth.

4.10 Installing a two-piece 400 mL reactor

- 1 Push the glass stirrer or the stirrer shaft with anchor or pitched-blade element through the central opening of the reactor cover before you place the cover onto the reactor.
- 2 Connect the stirrer shaft to the stirrer motor.
- 3 Place the reactor in the reactor holder.
- 4 Place the O-Ring on the cover rim and place the cover on the reactor.
- 5 Place the clamp chain carefully around the cover and the reactor flanges.
- 6 Screw the hook in or out so that the buckle can be easily closed (with one finger).

4.11 Turn on Device


- Power is connected.
- Cooling is connected and running.
- Purging of instrument is connected and running.
- Touchscreen is connected
- Press the ON/OFF button on the front side of the instrument.
 - ⇒ The touchscreen shows a splash screen during start up phase.
- ⇒ You can use the device as soon as the mainscreen appears.

5 Operation


All operations can be changed individually for reactor zone 1 or 2.

5.1 Select Reactor Type


1 Tap the button **Reactor 400 mL**.

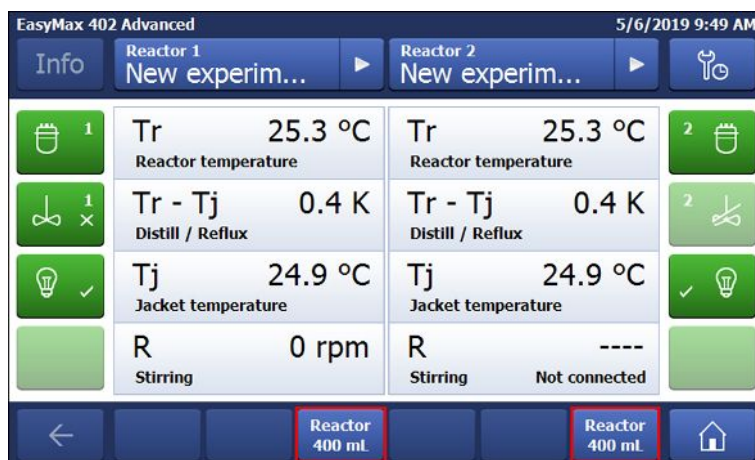

2 Tap the **Reactor type** field.

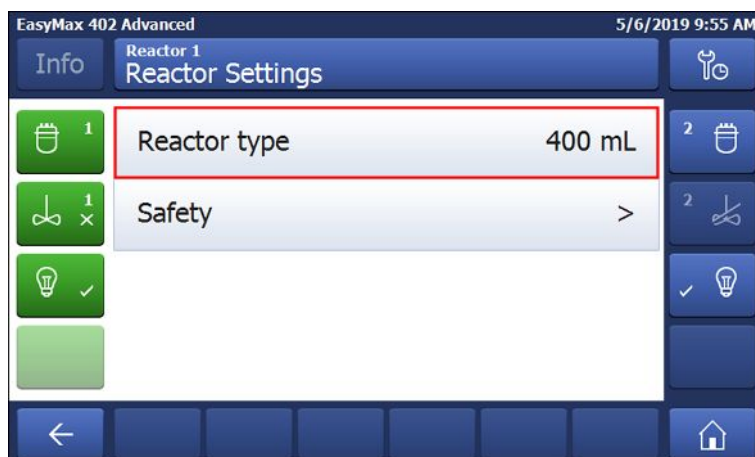



3 Select the reactor type that is installed.

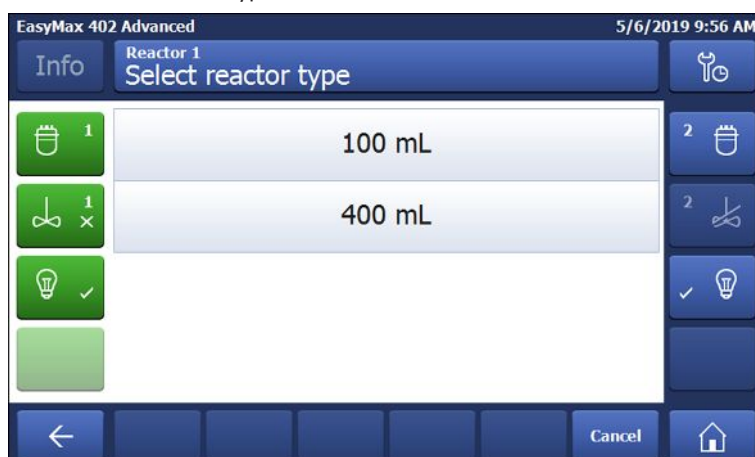

⇒ Make sure the safety settings for the reactor are still within the range.

5.2 Change Safety Settings


- 1 Tap the Reactor button.
- 2 Tap on the **Safety** field.

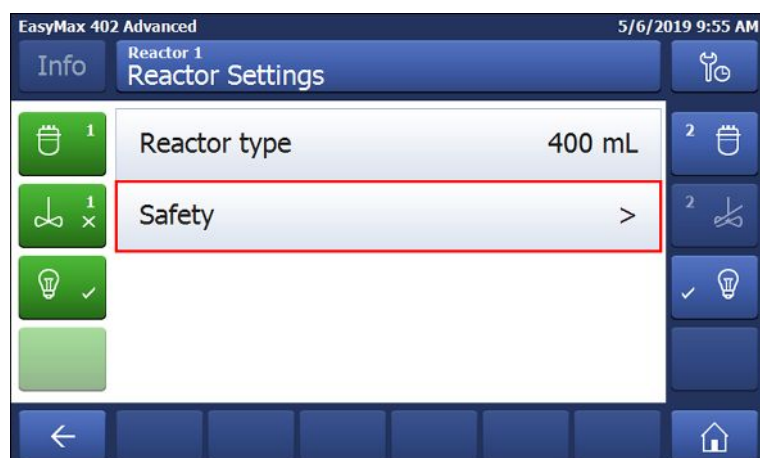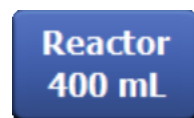


- 3 Change the necessary parameters according to your experiment and setup.

5.2.1 Change Safety Temperature (T safe)

- 1 Tap on **T safe**.

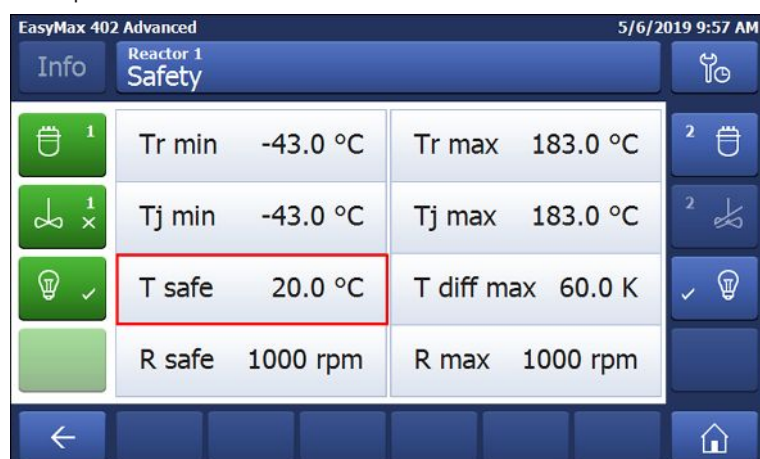

- 2 Enter a value for **T safe** that is valid for your experiment.

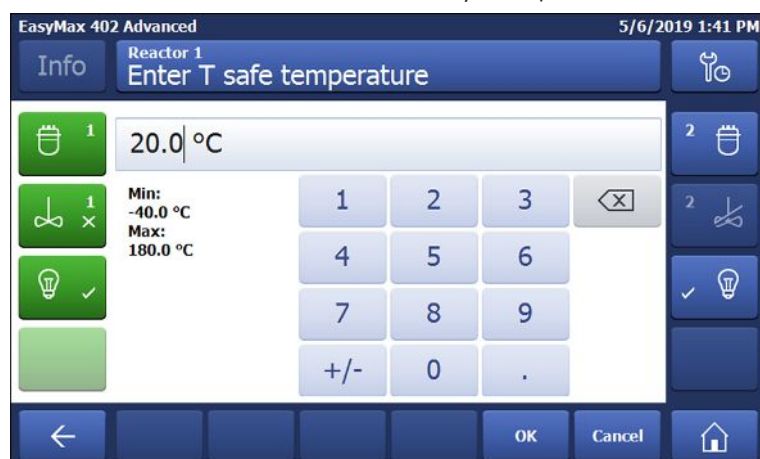

- 3 Tap **OK**.

5.2.2 Change Reaction Temperature limits (Tr)

- 1 Tap on **Tr max** or / and **Tr min**.

The screenshot shows the 'Reactor 1 Safety' screen of the EasyMax 402 Advanced control system. The interface is divided into a top header, a left sidebar with icons, and a main display area. The top header displays the system name and the date/time. The left sidebar contains icons for various functions: a calendar (1), a fan (1x), a lightbulb (checkmark), and an empty green box. The main display area is a table with two columns and four rows, showing setpoints for temperature and speed. The first two rows are highlighted with red borders. The right sidebar contains icons for a key (1), a fan (2), a scissors (2), a lightbulb (checkmark), and a house icon.

EasyMax 402 Advanced		5/6/2019 9:57 AM	
Reactor 1 Safety			
Tr min	-43.0 °C	Tr max	183.0 °C
Tj min	-43.0 °C	Tj max	183.0 °C
T safe	20.0 °C	T diff max	60.0 K
R safe	1000 rpm	R max	1000 rpm

- 2 Enter a value for **Tr max** and **Tr min** that is valid for your experiment.
- 3 Tap **OK**.

5.2.3 Change Range of Jacket Temperature (Tj)

- 1 Tap on **Tj min** or / and **Tj max**.

EasyMax 402 Advanced 5/6/2019 9:57 AM

Info Reactor 1 Safety

Tr min	-43.0 °C	Tr max	183.0 °C
Tj min	-43.0 °C	Tj max	183.0 °C
T safe	20.0 °C	T diff max	60.0 K
R safe	1000 rpm	R max	1000 rpm

- 2 Enter a value for **T_j min** and **T_j max** that is valid for your experiment.
- 3 Tap **OK**.

5.2.4 Change T diff max

- 1 Tap on **T diff max.**

EasyMax 402 Advanced

5/6/2019 9:57 AM

Info Reactor 1 Safety

1	Tr min	-43.0 °C	Tr max	183.0 °C	2
1 x	Tj min	-43.0 °C	Tj max	183.0 °C	2
✓	T safe	20.0 °C	T diff max	60.0 K	✓
	R safe	1000 rpm	R max	1000 rpm	

- 2 Enter a value for **T diff max** that is valid for your experiment.

- 3 Tap **OK**.

5.2.5 Change Rsafe

- 1 Tap on **Rsafe**.

- 2 Enter a value for **Rsafe** that is valid for your experiment.
- 3 Tap **OK**.


5.2.6 Change Rmax

- 1 Tap on **Rmax**.

- 2 Enter a value for **Rmax** that is valid for your experiments.
- 3 Tap **OK**.

5.3 Start an Experiment

- 1 Tap the experiment button on the main screen.

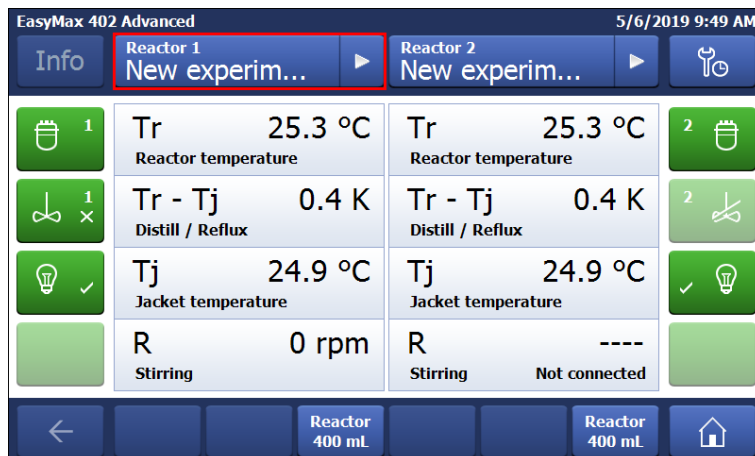


- 2 Enter an **experiment name**.
 - 3 Tap **Start** to start the experiment.
- ⇒ All tasks that are executed will be saved under the experiment and available for export.

5.4 Change Stirrer Speed

Note The value cannot be higher than the safety limit value.

- A stirrer is connected.
- The touchscreen is connected.

- 1 Tap on the **R** field.

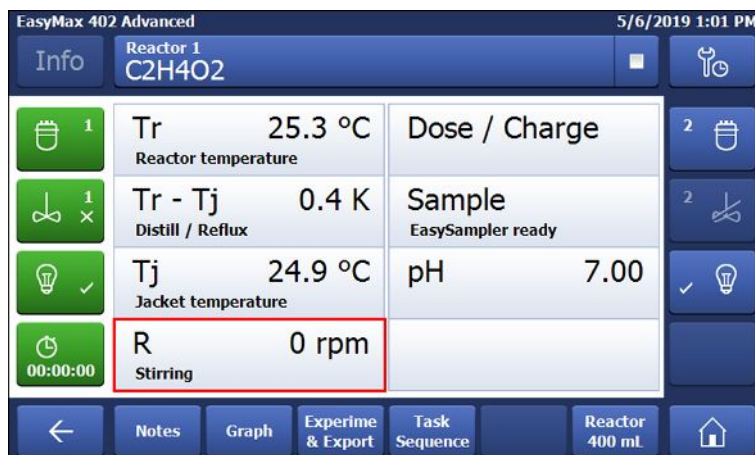


- 2 Enter the desired value.
 - 3 Tap **Start**.
- ⇒ The stirrer will immediately start stirring.

5.5 Change Tj

Note The value cannot be higher than the safety limit value.

- 1 Tap the **Tj** value field on the main screen.

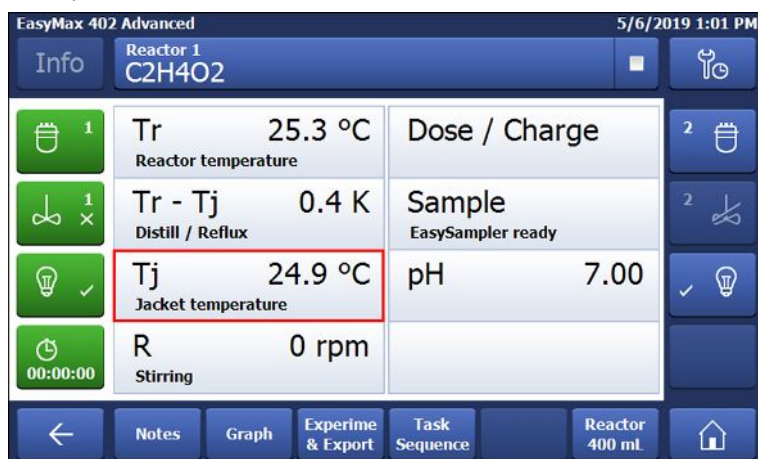


- 2 Enter the end temperature for **Tj**.
 - 3 Tap **Start** to initiate the task.
- ⇒ The task will start immediately.

5.6 Change Tr

Note The value cannot be higher than the safety limit value.

- A Tr sensor is connected to the thermostat.


- 1 Tap the **Tr** value field on the main screen.

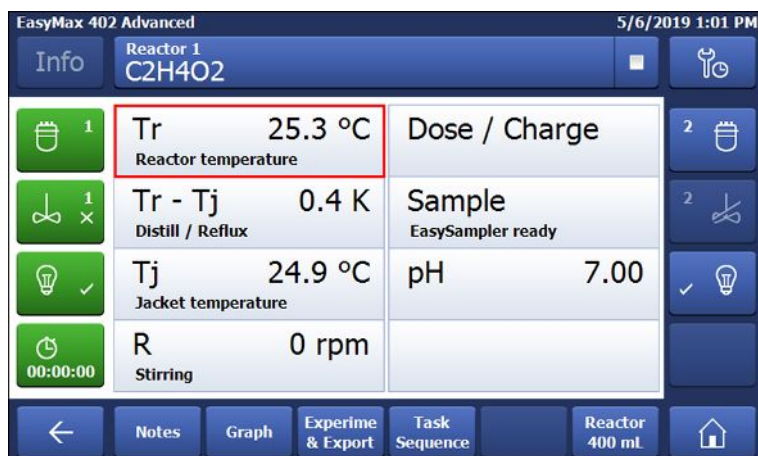

- 2 Enter the end temperature for Tr.
- 3 Tap **Start** to initiate the task.

5.7 End an Experiment

- 1 Tap on the Stop button on the mainscreen.

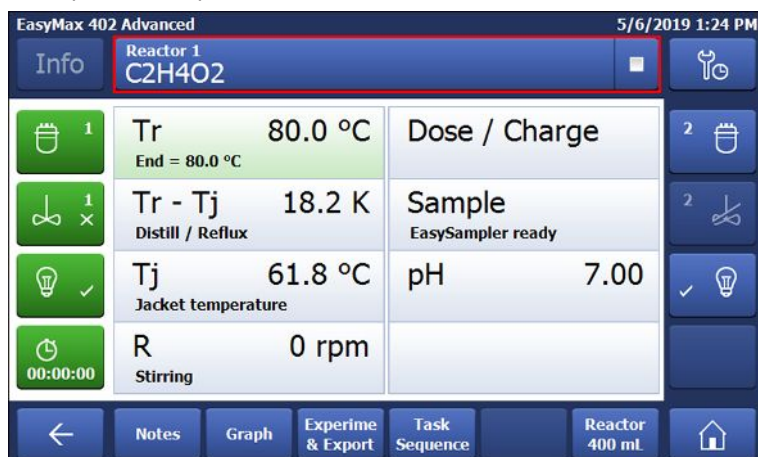

2 Select your preferred option for experiment end conditions.



3 Tap **OK**.

⇒ Your experiment is stored on the device and can be exported.

6 Maintenance

Maintenance tasks have to be performed in accordance with the instructions given in this chapter. After performing any maintenance tasks, it should be ensured that the device still meets all safety requirements.

6.1 Update Firmware

The latest firmware versions and instructions for installation are available on the following website:

<https://community.autochem.mt.com/?q=software>

6.2 Checking the Reactor

To check the reactor vessel for possible damage (scratches and cracks), it must be empty, clean, dry and open. Small hairline cracks can be detected by refraction using an additional light source (focused, not dispersed light).

6.3 Cleaning the Instrument



CAUTION

Hot instrument parts

Touching hot parts of the instrument can result in burnings.

- Do not clean the instrument before all parts have reached room temperature.



NOTICE

Damage to the device due to incompatible cleaning agents

Inappropriate cleaning agents could damage the housing of the device.

- 1 Use the described cleaning agent.
- 2 Should you use other cleaning agents, ensure that they are compatible with the housing material.

The housing of the instrument is not watertight (i.e. splash proof). We therefore recommend that you clean it with a damp cloth using ethanol.

If you have questions about the compatibility of cleaning agents, contact your authorized METTLER TOLEDO dealer or service representative.

6.4 Disposal

In conformance with the European Directive 2012/19/EU on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device. Should this device be passed on to other parties, the content of this regulation must also be related.



7 Technical Data

Certifications regarding this product can be found at <https://www.mt.com/us/en/home/search/compliance.html/>
The product name of your device is the model number.

Power supply

AC Power Adapter rating	Voltages	100 - 240 VAC
	Frequency	50 Hz / 60 Hz
	Permissible voltage fluctuation	±15%
Instrument rating	Power Consumption	Max. 1000 VA

Connections

All electrical connections	Not limited energy
USB	Support of USB 2.0
Electrical connectors	RS232, USB, CAN, Ethernet and touchscreen

Ambient Conditions

Humidity	Max. relative humidity 80 % for temperatures up to 31 °C decreasing linearly to 50 % relative humidity at 40 °C
Altitude	Up to 2000 m
Overvoltage category	II
Pollution degree	2
Ambient temperature	5 °C...40 °C
Usage	For indoor use only


Materials

Cover plate	Stainless steel coated with PFA/FEP
Housing material	Powder-Coated stainless steel
Connectors for purge gas	Stainless steel, nickel-plated brass
Purge gas lines	PVC, FEP, PP, PVDF, PTFE, aluminum
Coolant system (wetted parts)	PVC, PTFE, PVDF, copper
Cooling connector	Nickel-plated brass
Holder for lab bars	Aluminium
LEMO connectors for Tr sensor and overhead stirrer	Chrome-plated brass with protection cap in PSU
Reactor window	Borosilicate glass 3.3
Receptacles for reactors	Anodized aluminum
Fixing rings of thermostats	PTFE C25
USB connector	Stainless steel with protection cap in POM
On/Off switch	Stainless steel
Reactors	Borosilicate glass 3.3 and PTFE
Overhead stirrer shaft	Borosilicate glass 3.3, Alloy 22 or PTFE
Touchscreen	PA 12, aluminum
Protective cover for Touchscreen	Barex®

Device

Weight incl. Touchscreen	20 kg
Backlight	2 LED per reactor

Dimensions

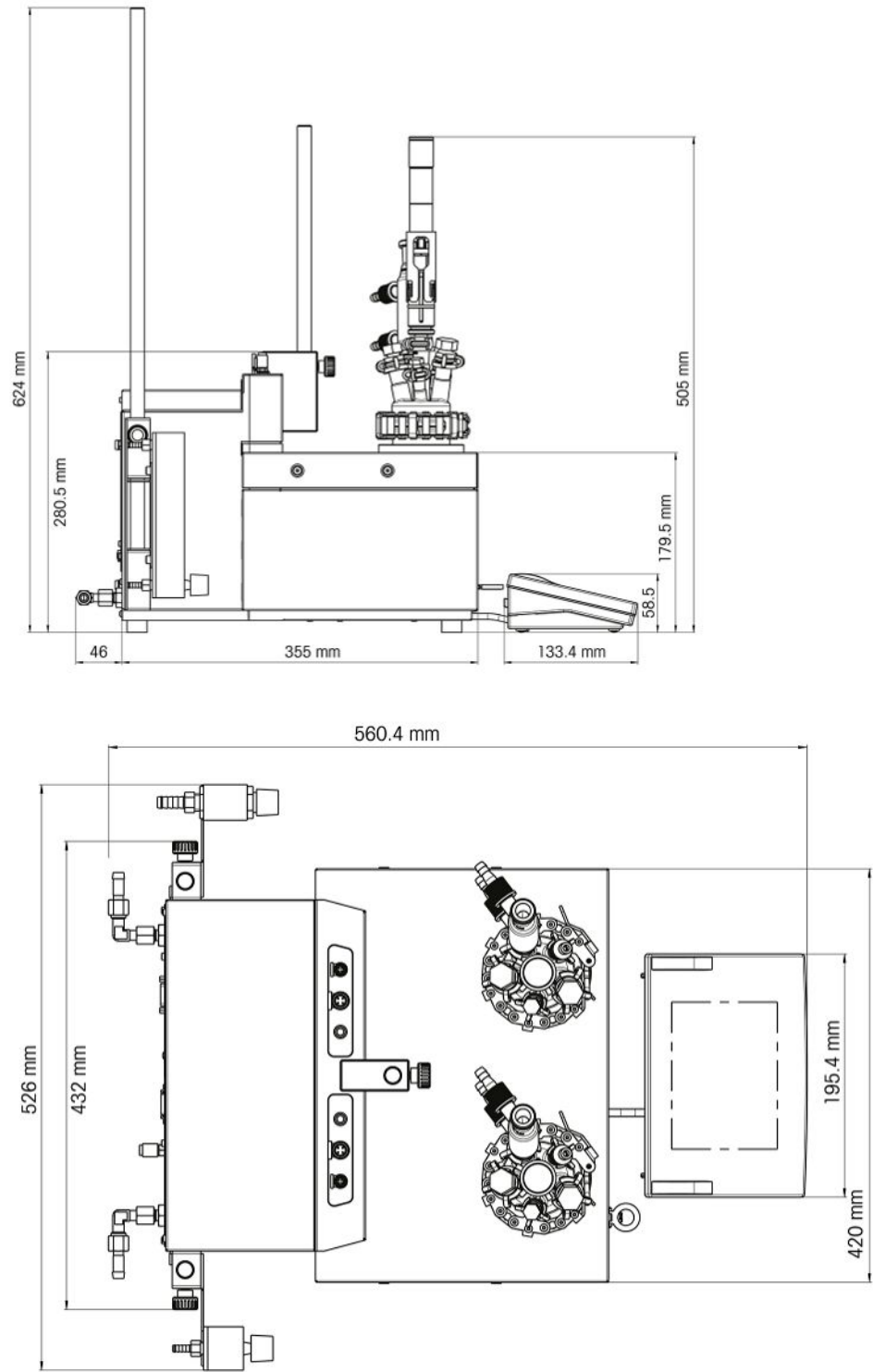

7.1 Thermostat

Power

Heating	360 W per thermostat
Cooling	Max. 150 W per thermostat

Temperatures

Range	Tj: - 40 °C (with Tc < -15 °C) to 180 °C Tr: - 40 °C (with Tc < -15 °C) to 180 °C* Tc: - 40 °C to 60 °C * The maximum and minimum Tr depends on heat transfer through the jacket and heat generated by the reaction.
Resolution	Tj: 0.1 K Tr: 0.1 K
Maximum permissible errors	1.0 °C for the whole range
Data recoding interval	Every 2 seconds

7.2 Reactors

Volumes	400 mL 100 mL
Pressure	0.05 bar to ambient pressure

Cover of reactors

100 mL glass cover	<ul style="list-style-type: none">• 2 x ST10/9• 3 x ST14/23<ul style="list-style-type: none">– 1x ST14/23: Central port for stirrer
100 mL PTFE cover	<ul style="list-style-type: none">• 1 x ST10/9• 3 x ST14/23<ul style="list-style-type: none">– 1x ST14/23: Central port for stirrer• 2 x ST19/26
100 mL PTFE 19 mm cover	<ul style="list-style-type: none">• 1 x ST10/19• 2 x ST14/23• 1 x 1/2" NPT• 1 x 19.25 mm• 1 x ST14 with 3/8" NPT: Central port for stirrer
400 mL glass cover	<ul style="list-style-type: none">• 4 x ST14/23<ul style="list-style-type: none">– 1x ST14/23: Central port for stirrer• 2 x ST19/26
400 mL PTFE cover	<ul style="list-style-type: none">• 3 x 1/2" NPT• 2 x 3/4" NPT• 1 x ST 14/35: Central port for stirrer

7.3 Stirrer

Overhead stirrer

Operating mode	Control to constant value or ramp
Data recording interval	Every 2 seconds
Speed range	50 to 1000 rpm

7.4 Cooling

Cooling medium	<ul style="list-style-type: none"> • Water (unpolluted); otherwise install a filter • Ethylene glycol • Ethanol • Silicone oil <p>Other cooling media have to be compatible with the wetted materials of the cooling system Technical Data</p>	
Min. flow of cooling media	2 L/min	
Cooling types	In-house coolant supply or cryostat	
Max. pressure of cooling media	<ul style="list-style-type: none"> • Without flowmeter: 3.5 bar • With flowmeter: 2 bar 	
Temperature (Tj)	In-house coolant supply	Cooling power of cryostat
> - 10 °C	At 15 °C	1000 W at 20 °C
- 40 °C	Not possible	Min. 450 W at -10 °C

7.5 Purge gas

Instrument purge gas

Max. inlet pressure	7 bar
Min. gas flow	As required by experiment

Reactor purge gas

Max. inlet pressure	7 bar
Min. gas flow	As required by experiment

To protect your product's future:

METTLER TOLEDO Service assures the quality, measuring accuracy and preservation of value of this product for years to come.

Please request full details about our attractive terms of service.

www.mt.com

For more information

Mettler-Toledo GmbH

Im Langacher 44
8606 Greifensee, Switzerland
www.mt.com/contact

Subject to technical changes.
© Mettler-Toledo GmbH 07/2019
30092611B



30092611